



Anton Paar

Measure
what is measurable
and make measurable
that which is not.

Galileo Galilei (1564-1642)

Instruction Manual and Safety Information

SurPASS 3
SurPASS 3 Eco

(Original Instruction)

Disclaimer

This document may contain errors and omissions. If you discover any such errors or if you would like to see more information in this document, please contact us at our address below. Anton Paar assumes no liability for any errors or omissions in this document.

Changes, copyright, trademarks etc.

This document and its contents may be changed or amended by Anton Paar at any time without prior notice.

All rights reserved (including translation). This document, or any part of it, may not be reproduced, changed, copied, or distributed by means of electronic systems in any form (print, photocopy, microfilm or any other process) without prior written permission by Anton Paar GmbH.

Trademarks, registered trademarks, trade names, etc. may be used in this document without being marked as such. They are the property of their respective owner.

Further information

Published and printed by Anton Paar GmbH, Austria

Copyright © 2023 Anton Paar GmbH, Graz, Austria

Address of the instrument producer:

Anton Paar GmbH
Anton-Paar-Str. 20
A-8054 Graz / Austria – Europe

Tel: +43 (0) 316 257-0

Fax: +43 (0) 316 257-257

E-Mail: info@anton-paar.com

Web: www.anton-paar.com

Date: January 2023

Document number: D85IB016EN-F

Instruction Manual and Safety Information

SurPASS 3 **SurPASS 3 Eco**

(Original Instruction)

Contents

1 Safety Instructions.....	5
1.1 General Safety Instructions.....	5
1.2 Conventions for Safety Messages	6
1.3 Safety Signs on the Instrument.....	6
2 SurPASS 3 - An Overview	8
2.1 Intended Use of the Instrument.....	8
2.2 Functional Elements	9
2.3 System Requirements.....	10
2.4 Water Quality and Measurement Conditions	10
3 Preparing the Instrument	12
3.1 Placing the SurPASS 3.....	12
3.2 Removing the Transport Lock.....	12
3.3 Mains Connection	13
3.4 PC Connection.....	14
3.5 Mounting the Electrolyte Hoses	14
3.6 Connecting the Conductivity and pH Electrodes.....	14
3.7 Mounting the Hoses of the Titration Unit (not applicable to SurPASS 3 Eco).....	15
3.8 SurPASS 3 Accessory Kit for Liquid Handling	15
3.9 SurPASS 3 Pressure Control Unit (not applicable to SurPASS 3 Eco)	16
3.10 SurPASS 3 Software.....	17
4 Putting SurPASS 3 into Operation	20
4.1 pH Electrode Calibration	20
4.2 Conductivity Electrode Calibration	21
4.3 Filling of the Titration Unit (not applicable to SurPASS 3 Eco)	22
4.4 SurPASS 3 Verification Measurement with Cotton Fabrics	22
5 Performing a Measurement.....	27
5.1 Mounting the Measuring Cells on SurPASS 3	27
5.2 Rinsing the System	28
5.3 Starting a Zeta Potential Measurement	29
5.4 Adsorption Kinetics Measurement (not applicable to SurPASS 3 Eco)	31
5.5 End of Measurement.....	31
5.6 Data Analysis	33
6 Upkeep and Cleaning	35
6.1 Software Administration	35
6.2 Upkeep and Cleaning after Standard Operation.....	35
6.3 Upkeep and Cleaning after Adsorption Operation	35
6.4 Cleaning and Drying the Measuring Cells.....	36
6.5 Wetted Parts	36
6.6 Exchange of the Electrodes	36
6.7 Manual Cleaning of the Electrolyte Circuit.....	36
6.8 Removing and Mounting of the Syringes of the Titration Unit (not applicable to SurPASS 3 Eco).....	37
6.9 Venting the Gas Supply Line (not applicable to SurPASS 3 Eco)	37
6.10 Storing and Transporting the Instrument	37
6.11 Packing the Instrument for Returns	37
7 Maintenance and Repair.....	38
7.1 Maintenance Performed by an Authorized Anton Paar Service Engineer	38
7.2 Repair Performed by an Authorized Anton Paar Representative	38
Appendix A: Technical Data	39
Appendix B: Declaration of Conformity.....	41
Appendix C: Requirements for Nitrogen Supply	42

1 Safety Instructions

- Read the documentation before using the SurPASS 3 instrument.
- Follow all hints and instructions contained in the documentation to ensure the correct use and safe functioning of the SurPASS 3 instrument.
- The documentation is a part of the product. Keep this document for the complete working life of the product and make sure it is easily accessible to all people involved with the product. If you receive any additions or revisions to the documentation from Anton Paar GmbH, these must be treated as part of the documentation.
- Continue operating SurPASS 3. Do not operate the instrument under conditions which could result in damage to goods and/or injuries and loss of life.
- If you observe leakage underneath the SurPASS 3 instrument, immediately stop the measurement and disconnect the SurPASS 3 instrument from its power supply. Identify the reason for leakage and contact your Anton Paar representative.
- Check SurPASS 3 for chemical resistance to the samples and cleaning agents. See section 6.5 Wetted Parts for a list of materials that come in contact with the electrolyte solution and titration liquids.

1.1 General Safety Instructions

Liability

- The documentation does not claim to address all safety issues associated with the use of the instrument and samples. It is your responsibility to establish health and safety practices and determine the applicability of regulatory limitations.
- Anton Paar GmbH only warrants the proper functioning of SurPASS 3 if no adjustments have been made to the mechanics, electronics, and firmware. Do not perform any modifications on the SurPASS 3 instrument.
- Only use SurPASS 3 for the purpose described in the documentation. Anton Paar GmbH is not liable for damages caused by incorrect use of SurPASS 3.

Installation and use

- SurPASS 3 is **not** an explosion-proof instrument and therefore must not be operated in areas with risk of explosion.
- The installation procedure should only be carried out by authorized personnel who are familiar with the installation instructions.
- Do not use any accessories or spare parts other than those supplied or approved by Anton Paar GmbH.
- Make sure all operators are trained to use the instrument safely and correctly before starting any applicable operations.
- In case of damage or malfunction, do not con-

Maintenance and service

- The results delivered by SurPASS 3 not only depend on the correct functioning of the instrument, but also on various other factors. We therefore recommend you have the results checked (e.g. plausibility tested) by skilled personnel before consequential actions are taken based on the results.
- Service and repair procedures may only be carried out by authorized personnel or by Anton Paar GmbH.

Disposal

- Concerning the disposal of SurPASS 3, observe the legal requirements in your country.

Returns

- For repairs send the cleaned SurPASS 3 to your Anton Paar representative. Only return the instrument together with the filled out RMA (Return Material Authorization) and the form "Safety Declaration for Instrument Repairs". Please download the Safety Declaration form from our website www.anton-paar.com.
- Do not return instruments which are contaminated by radioactive materials, infectious agents or other harmful substances that cause health hazards.

Precautions for highly inflammable samples and cleaning agents

- Observe and adhere to your national safety regulations for handling the measured samples

1 Safety Instructions

(e.g. use of safety goggles, gloves, respiratory protection etc.).

- Only store the minimum required amount of sample, cleaning agents and other inflammable materials near SurPASS 3.
- Do not use inflammable, toxic or explosive samples with SurPASS 3.
- Do not spill sample/cleaning agents or leave their containers uncovered. Immediately remove spilled sample/cleaning agents.
- Make sure that the setup location is sufficiently ventilated. The environment of SurPASS 3 must be kept free of inflammable gases and vapors.
- Connect SurPASS 3 to the mains via a safety switch located at a safe distance from the instrument. In an emergency, turn off the power using this switch instead of the power switch on SurPASS 3.
- Supply a fire extinguisher.
- Ensure the sufficient supervision of SurPASS 3 during operation.

1.2 Conventions for Safety Messages

The following conventions for safety messages are used in this instruction manual:



WARNING

Description of risk.

Warning indicates a hazardous situation which, if not avoided, could result in death or serious injury.



CAUTION

Description of risk.

Caution indicates a hazardous situation which, if not avoided, could result in minor or moderate injury.

NOTICE

Description of risk.

Notice indicates a situation which, if not avoided, could result in damage to property.



CAUTION

Electrostatic sensitive device

Caution indicates a situation which, if not avoided, could result in damage to property.



Wear protective gloves

Wear protective gloves when handling the SurPASS 3 instrument.



Wear safety goggles

Wear safety goggles when handling the SurPASS 3 instrument.



Use protective clothing

Use protective clothing when handling the SurPASS 3 instrument.

TIP: Tip gives extra information about the situation at hand.

1.3 Safety Signs on the Instrument

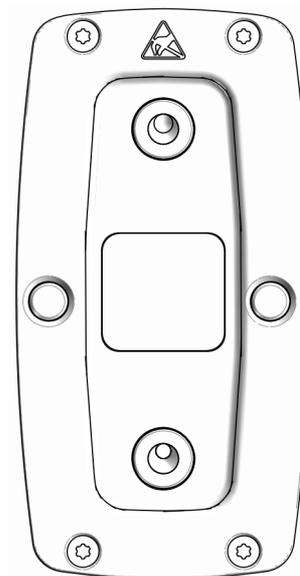


Fig. 1: Warning signs on the instrument where the measuring cells are to be connected



WARNING

Hazardous liquids

The electrolyte circuit may contain hazardous liquids. Make sure to empty the electrolyte circuit before removing the measuring cell.

**CAUTION*****Electrostatic sensitive device***

SurPASS 3 is an ESD sensitive device. Do not touch the electrodes.

**WARNING*****Hazardous liquids***

The titration unit of SurPASS 3 is intended for use with acid and base. Always empty the titration unit and rinse it before removing the titration hoses.

NOTICE

Take care that the warning symbols remain clearly legible.

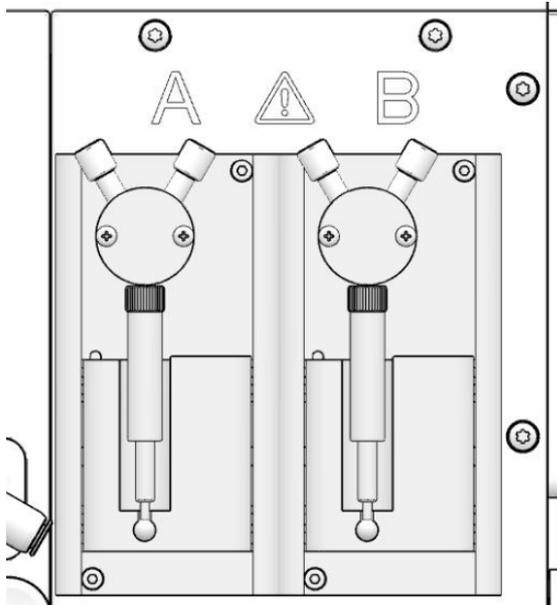


Fig. 2: Warning sign on the SurPASS 3 titration unit (not applicable to SurPASS 3 Eco)

2 SurPASS 3 - An Overview

Thank you for buying SurPASS 3 for the zeta potential analysis of **macroscopic** solid surfaces.

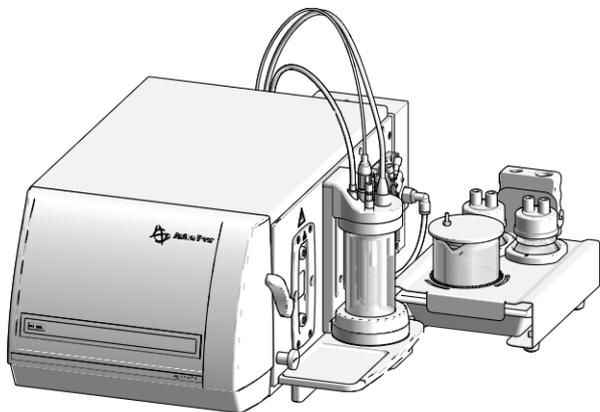


Fig. 3: SurPASS 3 instrument

The SurPASS 3 instrument is available in two different configurations: As fully equipped SurPASS 3 for high-end zeta potential analysis as it is described in this instruction manual, and as entry-level device for single-point measurements, SurPASS 3 Eco. With SurPASS 3 Eco the recording of automated pH dependences and studies of adsorption kinetics are disabled. Please note that SurPASS 3 Eco is upgradeable to a fully equipped SurPASS 3 on-site by an Anton Paar certified service engineer, in case you wish to exploit the full potential of your SurPASS 3 instrument in the near future.

TIP: In this instruction manual, instrument and software features which are not accessible with SurPASS 3 Eco, are labeled as such.

Together with the SurPASS 3 instrument you have received the SurPASS 3 software for the remote control of SurPASS 3 with a PC or laptop computer.

The SurPASS 3 software offers the following functions:

- **Remote control of the SurPASS 3 instrument**
The used measuring cell is automatically detected and the zeta potential analysis can be started at the click of a button. All key parameters are displayed in a single view.

- **Automatic detection of the isoelectric point** (*not applicable to SurPASS 3 Eco*)
SurPASS 3 detects whether acidic or alkaline titration has to be applied for the automated determination of the isoelectric point of a sample.
- **Automatic pH scans** (*not applicable to SurPASS 3 Eco*)
Complete pH titrations can be performed fully automatically over a desired pH range.
- **Adsorption kinetics** (*not applicable to SurPASS 3 Eco*)
Time-resolved adsorption studies from the liquid phase onto solid surfaces can be performed by monitoring the change in streaming potential or streaming current while changing the composition of the electrolyte solution.
- **Display and export of data**
Measurement data can be loaded and displayed at any time. It is also possible to compare measurement results of different files. Furthermore, all data are automatically saved for user-specific evaluation in a spreadsheet software.

2.1 Intended Use of the Instrument

The SurPASS 3 instrument is applicable for the zeta potential analysis of solids of almost any shape and size like fibers, foils, sheets, textiles, powders or granules. The principle for the zeta potential analysis with the SurPASS 3 instrument is based on the streaming potential and streaming current method.

In combination with the integrated titration unit, it is possible to investigate the dependence of the zeta potential on the pH fully automatically, thus determining the isoelectric point of a sample surface.

The SurPASS 3 instrument is intended for use with dilute aqueous solutions of inorganic salts, acids and bases. For standard measurements, the background electrolyte solution composes of a 0.001 mol/l aqueous NaCl or KCl solution. For adjusting the pH value of the background electrolyte solution, the usage of 0.05 mol/l HCl and 0.05 mol/l NaOH or KOH, respectively, is recommended.

2.2 Functional Elements

2.2.1 Front View

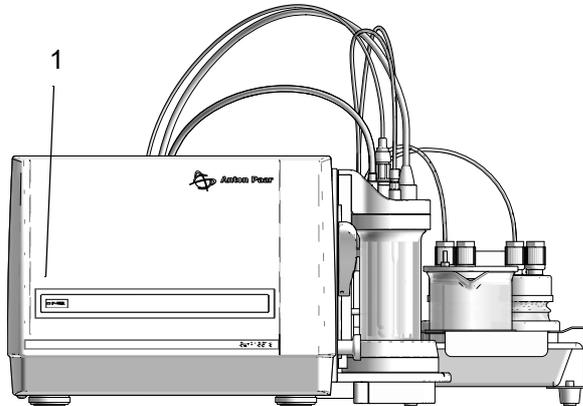


Fig. 4: SurPASS 3 front view. The green LED (1) indicates that the instrument is switched on.

2.2.2 Rear View

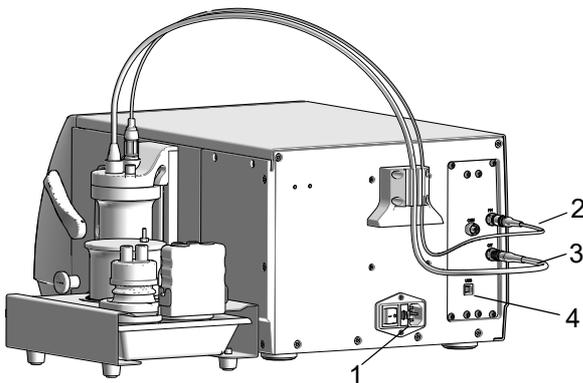


Fig. 5: SurPASS 3 Eco rear view

- 1 power switch
- 2 connector for pH electrode
- 3 connector for conductivity electrode
- 4 connector for USB cable

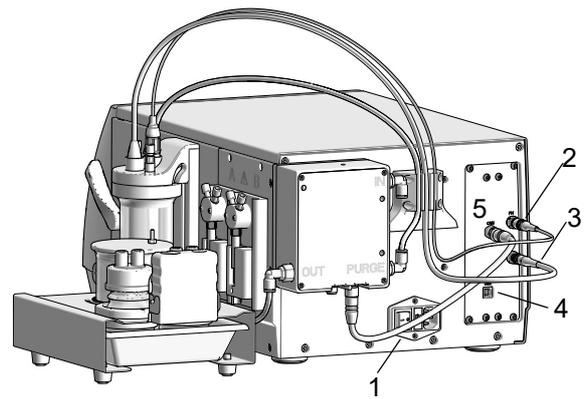


Fig. 6: SurPASS 3 rear view

- 1 power switch
- 2 connector for pH electrode
- 3 connector for conductivity electrode
- 4 connector for USB cable
- 5 connector for CAN cable for pressure control unit

2.2.3 Right Side View

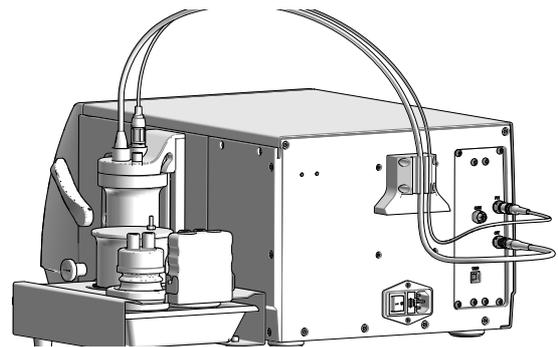


Fig. 7: Right side view of SurPASS 3 Eco

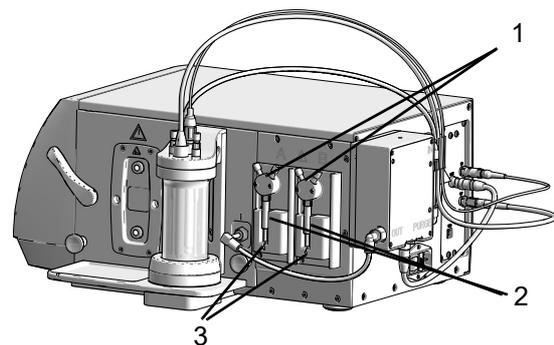


Fig. 8: Right side view of SurPASS 3

- 1 3-way valve of SurPASS 3 titration unit
- 2 1.0 ml syringe of SurPASS 3 titration unit
- 3 piston sled of SurPASS 3 titration unit

2.3 System Requirements

The following hardware configuration is required to successfully install and run the SurPASS 3 software on your computer:

- Intel® Core™ i3 CPU
- 4 GB RAM (depends on the operating system)
- Screen resolution of at least 1024 x 768 at True Color (32 bit), best experience with a monitor with an aspect ratio of 16:9 or 16:10
- Use of a dedicated graphics card that supports Direct X 9 is recommended
- 20 GB free disk space on system drive
- Mouse and keyboard

The SurPASS 3 software can be installed on one of the following operating systems:

- Microsoft Windows® 7
 - 32 bit and 64 bit are supported
 - Service Pack 1 or later is required
 - Windows 7 Starter Edition is not supported
- Microsoft Windows® 8
 - 32 bit and 64 bit are supported

- Microsoft Windows® **Server 2008 / 2008 R2**
 - 32 bit and 64 bit are supported
 - Service Pack 2 or higher is required for Windows Server 2008
- Microsoft Windows® **10**
 - 32 bit and 64 bit are supported

NOTICE

Please make sure that all requirements are fulfilled to guarantee trouble-free operation of the SurPASS 3 software.

2.4 Water Quality and Measurement Conditions

The zeta potential is very sensitive to the outermost solid surface. Care needs to be taken when it comes to the selection of the chemicals used for measurement. High purity chemicals need to be used for measurement, since contamination may affect the analyses especially of hydrophobic samples or samples with a small surface area (e.g. polymer films).

Table 1: Required chemicals for successful SurPASS 3 measurements

ultrapure water (ASTM I)	for rinsing the electrolyte hoses and for preparing the electrolyte solution for measurement
KCl dry salt, > 99.5 % <i>p.a.</i> or NaCl dry salt, > 99.5 % <i>p.a.</i>	for the preparation of 0.001 mol/l electrolyte solution
0.05 mol/l HCl*	acid for pH titration
0.05 mol/l NaOH or KOH*	base for pH titration
three buffer solutions (e.g. pH 3, 7, 10)	for pH calibration
0.1 mol/l KCl solution or conductivity standard	for conductivity calibration
* A concentration of 0.05 mol/l is best suited for the pH titration in the range of pH 3-9. If the zeta potential analysis is required at lower or higher pH, the concentration of the stock solution of acid and base must be increased accordingly.	

A crucial requirement for a successful measurement is the quality of water, which is used for the operation of the SurPASS 3 instrument. The water used must comply with the American Standard for Testing and Materials I (ASTM I).

Table 2: Minimum ASTM I standard

conductivity (at 25 °C)	0.056 $\mu\text{S}/\text{cm}$
electric resistivity (at 25 °C)	18.0 $\text{M}\Omega \text{ cm}$
TOC	max. 100 $\mu\text{g}/\text{l}$
pH value*	5.5 - 6.5
* not specified by ASTM I	

NOTICE

- It is not possible to perform successful SurPASS 3 verification measurements (see chapter 4.4) if the water quality is not sufficient. In the worst case, the water used for measurement can even seriously contaminate the electrolyte circuit.
- Commercial water purification units are specified in terms of ASTM standards but only provide the indicated water quality if they are also maintained regularly, in intervals as recommended by the manufacturer.
- The use of bottled ultrapure water is possible if no appropriate water purification unit is available.
- It is recommended to continuously purge the electrolyte solution with nitrogen 5.0 throughout the duration of the measurements. Do not use compressed air or inert gas with a lower quality.

NaCl or KCl used for the fresh preparation of the measuring electrolyte needs to be *pro analysis* grade. It is important that the salt is dry and not contaminated by any other chemical. It is recommended to use fresh NaCl *p.a.* or KCl *p.a.* for preparing the measuring electrolyte solution by putting 15 mg of KCl salt (12 mg of NaCl) and dissolving them in 200 ml ultrapure water. The preparation of stock solutions is not recommended. It is not recommended to use ready-made NaCl or KCl solutions, as they may contain stabilizers which also affect zeta potential measurements.

Acid and base come in contact with the electrolyte system and the sample during pH titration. It is recommended to exchange the 0.05 mol/l HCl and 0.05 mol/l NaOH (or KOH) solutions after every month.



WARNING

Risk of injury

Acids and bases may only be handled by skilled personnel who are aware of all risks associated to the used liquids. Refer to the reference guide D85IB016EN for the safety data sheets of 0.1 mol/l HCl and 0.1 mol/l NaOH.



Wear protective gloves

Wear protective gloves when handling acids and bases.



Wear safety goggles

Wear safety goggles when handling acids and bases.



Use protective clothing

Use protective clothing when handling acids and bases.

3 Preparing the Instrument

3.1 Placing the SurPASS 3

SurPASS 3 is a benchtop instrument. Select a suitable place for your SurPASS 3 instrument to avoid incorrect measurements.

- Make sure that the SurPASS 3 is placed on a table in a horizontal position to avoid an overflow of electrolyte from the collecting dike in case of malfunction.
- Make sure that the SurPASS 3 is placed in a way to reach the power switch easily.
- Do not place SurPASS 3 near heat sources.
- Avoid direct exposure of SurPASS 3 to sunlight.
- Prevent influence by vibrations.

TIP: When moving SurPASS 3, lift the instrument at the front and rear bottom.

3.2 Removing the Transport Lock

NOTICE

Before putting the SurPASS 3 instrument in operation, you need to connect the cable of the hydraulic valve motor and remove the transport lock (Kapton foil) that protects the hydraulic valve.

If you fail to do so before starting a process (e.g. Rinse), the SurPASS 3 software will display an error message:

Rinse failed

The operation failed on SurPASS 3
Error code: 02100004h

Please contact customer support and provide the error details.

Operation type: 4 'Current' sequence
Operation error code: -1 'Not translated!'
Operation status code: 8 'Sequence failed'

Instructions for removing the transport lock:

1. Make sure the SurPASS 3 instrument is

switched off.

2. Remove the two knurled screws in the corners of the front plate.



Fig. 9: Front plate of the plug-in module

3. Pull out the plug-in module completely.

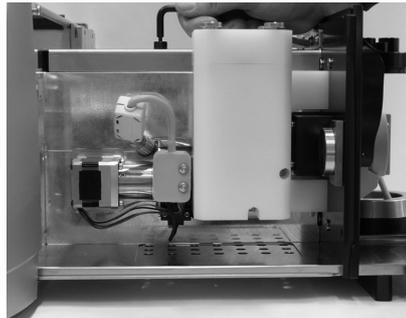


Fig. 10: Plug-in module

4. Take the plug of the hydraulic valve motor from its support in the Kapton foil.

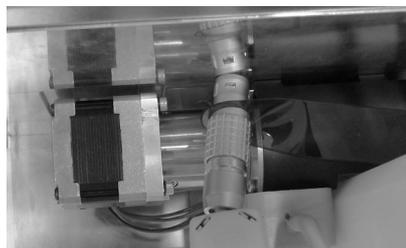


Fig. 11: Removing the Kapton foil from the plug

5. Connect the plug of the hydraulic valve motor to the connector on the back housing.



Fig. 12: Connecting the hydraulic valve motor

6. Insert the blade of the T20 Torx T-handle screw driver into the corresponding holes on the back side of the plug-in module housing. Support the valve motor and open the two screws to a small extent, such that the Kapton foil gets movable.



Fig. 13: Loosen the screws of the hydraulic valve motor

7. Pull out the Kapton foil that protects the hydraulic valve. Make sure the Kapton foil is pulled out as a whole. Fix the two screws again, hand-tight only.



Fig. 14: Pulling out the Kapton foil

8. Insert the plug-in module into the SurPASS 3 housing.
9. Lock the front panel by fixing the knurled screws on the bottom of the front panel.

NOTICE

The Kapton foil must NOT be discarded. If the instrument is kept out of operation for an extended period, the Kapton foil must again be placed between the ceramic discs - refer to section 9.7.4 in the reference guide D85IB016EN.

3.3 Mains Connection

The power cable receptacle and the power switch are located at the rear of the SurPASS 3 instrument (pos. 1 in Fig. 6). The mains voltage is indicated at the back side of SurPASS 3.



CAUTION

Mains Voltage

SurPASS 3 is designed for operation on mains fulfilling the following requirements:
115/230 VAC, 50/60 Hz, $\pm 10\%$ allowable mains voltage tolerance.



WARNING

Risk of injury

If you notice that the power cable is defective, immediately disconnect SurPASS 3 from the power supply. Replace the power cable.

3.3.1 Protective Grounding

Grounding of the SurPASS 3 instrument is done through a three-wire power cable which has to be connected to a protective ground connector socket.

No other kind of grounding is permitted.



WARNING

Risk of injury

When moving SurPASS 3 from cold to warm surroundings, parts may become conductive due to water condensation. Strictly obey all grounding regulations!

Interrupting the protective grounding inside or outside of SurPASS 3 is inevitably dangerous. Never use extension cables without a protective ground connector.

3.4 PC Connection

The SurPASS 3 instrument is controlled by a PC with the Microsoft Windows® based SurPASS 3 software. SurPASS 3 communicates with the PC via a USB cable which is connected at the rear side of SurPASS 3 in the socket *USB* (pos. 4 in Fig. 6).

3.5 Mounting the Electrolyte Hoses

The electrolyte hoses inside the SurPASS 3 instrument have already been mounted before shipment.

1. Pull the loose ends of the two hoses through the corresponding holes of the white beaker cover.

NOTICE

The beaker cover has been designed for use with a 250 ml glass beaker (60 mm diameter) which is supplied with the SurPASS 3 instrument.

3.6 Connecting the Conductivity and pH Electrodes

The conductivity electrode is equipped with a tem-

perature sensor and a single connector cable.

1. Plug the conductivity electrode cable in the socket *C/T* of the connector panel on the backside of the SurPASS 3 instrument (pos. 3 in Fig. 6).
2. Place the conductivity electrode in an appropriate hole of the beaker cover (pos. 2 in Fig. 15).



Fig. 15: Drawing of the beaker cover

- 1 Position for pH probe
- 2 Position for conductivity electrode
- 3 Positions (for acid and base dispense) for titration hoses
- 4 Position for nitrogen purge hose with throttle valve and electrolyte inlet

Proceed accordingly with the pH electrode:

1. Plug the pH electrode cable in the socket *PH* of the connector panel on the backside of SurPASS 3 (pos. 1 in Fig. 6).
2. Place the pH electrode in an appropriate hole of the beaker cover (pos. 1 in Fig. 15).

3.7 Mounting the Hoses of the Titration Unit (not applicable to SurPASS 3 Eco)

Two syringe pumps (syringe pump A and B) for automated dispense of titration liquid are integrated in the SurPASS 3 instrument. The corresponding 1 ml syringes are already mounted on the SurPASS 3 upon delivery. Follow the instructions given in chapter 6.8 if you need to remove and remount them for e.g. maintenance purposes or replacement.

Two sets of hoses are supplied with the SurPASS 3 accessories:

- 2 dispense lines (red and blue) with connector and ferrules on one end, and burette tip on the other end,
- 2 uptake lines (red and blue) with connector and ferrules on one end only



WARNING

Risk of injury

Handle the hoses for the titration unit only when they are clean and empty.

These hoses need to be mounted on the top left and right positions of the 3-way valves of the integrated titration unit.

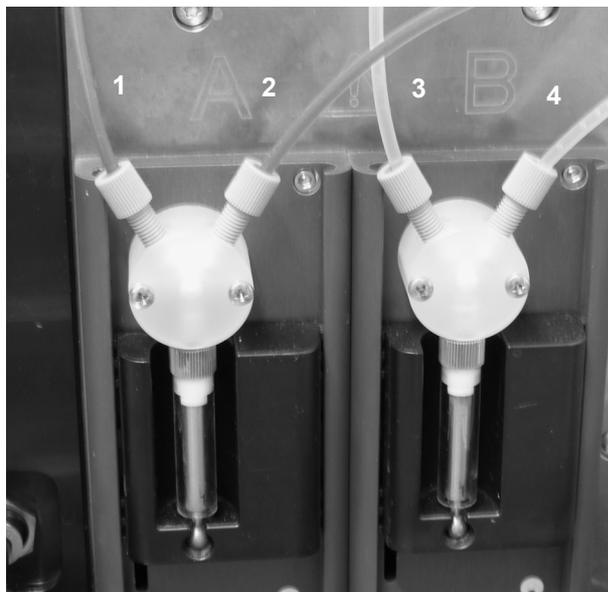


Fig. 16: Mounted hoses of the titration unit

- 1 Dispense line red with burette tip
- 2 Uptake line red (open end)
- 3 Dispense line blue with burette tip
- 4 Uptake line blue (open end)

1. Mount the nut of the dispense line in the top-left hole with thread of a 3-way valve.
2. Place the burette tip of a dispense line in the appropriate hole of the beaker cover (pos. 3 in Fig. 15)
3. Mount the nut of an uptake line in the top right hole with thread of a 3-way valve.
4. Fix the open end of the uptake line in the supply bottle with the titration liquid.



WARNING

Risk of injury

The SurPASS 3 instrument is intended for use with dilute aqueous solutions of inorganic salts, acids and bases. The usage of 0.05 mol/l HCl and 0.05 mol/l NaOH, respectively, is recommended for adjusting the pH value of the background electrolyte solution.

Always wear protective clothing, protective gloves and safety goggles when handling acids and bases.

TIP: To use the settings in the SurPASS 3 software, connect acid to syringe pump A and base to pump B.

5. Proceed accordingly with the second titration unit.



CAUTION

Hose connections

For a safe operation check the titration hoses and connections regularly.

3.8 SurPASS 3 Accessory Kit for Liquid Handling

The SurPASS 3 accessory kit for liquid handling provides a conscientious and safe supply of the used liquids. This accessory arranges the two supply bottles for the titration liquids, a 400 ml waste beaker with cover for connecting the outlet hose and two storage containers for the pH and conductivity electrodes, respectively.

Mount the SurPASS 3 accessory kit for liquid handling as shown in Fig. 17. Place the device on the right side of the SurPASS 3 instrument close to the titration unit. Make sure the hoses of the titration unit are mounted as described in chapter 3.7.

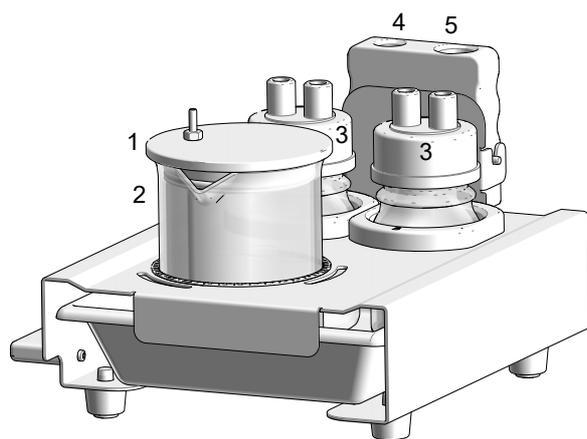


Fig. 17: SurPASS 3 accessory kit for liquid handling

- 1 Cover of waste beaker
- 2 Waste beaker
- 3 Storage bottle for acid, base
- 4 Storage position for the conductivity electrode
- 5 Storage position for the pH electrode



WARNING

Risk of injury

To avoid a flooding after the clean cycle, empty the waste beaker and the dike if there is any liquid in it.

3.9 SurPASS 3 Pressure Control Unit (not applicable to SurPASS 3 Eco)



WARNING

Risk of injury

- Always switch off the instrument before handling the hoses of the gas connection.
- After activation of the gas connection, the gas hose which connects the SurPASS 3 instrument with your in-house gas supply will remain pressurized. Make sure that all hoses are unpressurized before disconnecting any of the hoses (see chapter 6.9).

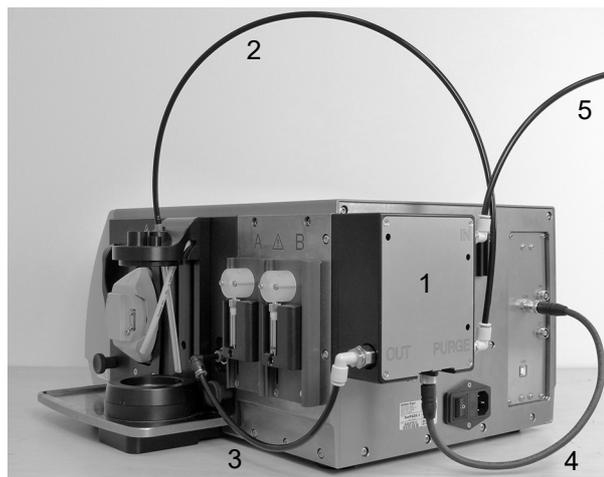


Fig. 18: SurPASS 3 external pressure control

- 1 Pressure control unit
- 2 Gas hose for purging the electrolyte solution
- 3 Gas hose for internal pressure supply
- 4 CAN cable for pressure control unit
- 5 Gas hose for nitrogen supply

1. Connect the end of hose (2) with the orange throttle valve (in direction of the arrow towards the beaker) to the electrolyte inlet to the T-connector. The other end of hose (2) still needs to be connected to the position "PURGE" of the pressure control unit (1). Mount the hose tightly into the connector by pushing it in until the end stop.
2. Upon delivery of the SurPASS 3 instrument, hose (3) is already connected between the position "OUT" of the pressure control unit (1) and the gas hose connector (position "p") on the instrument.
3. The CAN cable (4) connects the right one of the two sockets at the lower position of the pressure control unit (1) with the socket "CAN" of the connector panel on the backside of the SurPASS 3 instrument (pos. (5) in Fig. 6).
4. Hose (5) needs to be connected between the position "IN" on the pressure control unit (1) and the nitrogen supply in your laboratory. Refer to appendix C for details on the requirements for nitrogen supply and gas consumption.

TIP: Make sure that the nitrogen bubbles gently during the measurements to avoid the dissolution of the CO₂ from air.

**WARNING****Risk of injury**

As there is a risk of suffocation when using nitrogen for measurement and to purge the electrolyte solution, ensure adequate ventilation of the room.

NOTICE

The operating pressure of the pressure control unit, *i.e.* the pressure of the nitrogen supply in your laboratory, shall be set to 6 bar. The operating pressure must not exceed 10 bar. The outgoing pressure is adjusted in the software.

3.10 SurPASS 3 Software

3.10.1 Installation of the SurPASS 3 Software

For the installation of the SurPASS 3 software start the installation file and follow the instructions of the setup program. Before installing the software, read the system requirements and make sure that your PC or laptop computer meets the hardware and software requirements (see chapter 2.3).

TIP: *You must have administrator privileges for the installation of the SurPASS 3 software.*

NOTICE

Do not install this product on servers or workstations that run mission-critical applications.

The following software components will be automatically installed upon first installation of the SurPASS 3 software:

- Microsoft Windows Installer 4.5
- Microsoft NET Framework 4.5.1

You do not have to enter a license key for operation of SurPASS 3 with the SurPASS 3 software.

3.10.2 Graphical User Interface

Open the SurPASS 3 software by double-clicking on the icon, which was created on the PC desktop during installation.

The SurPASS 3 software is a Microsoft Windows®

based software offering the standard features of Windows programs.

Software functions specific for the operation of the SurPASS 3 instrument are (numbers in brackets refer to Fig. 19) described below:

TIP: *Certain software features are disabled if you are operating a SurPASS 3 Eco.*

- **Application Button (1):** Click on the application button <Sp> to open the *settings view*.
- **Settings View:** In the *settings view* you can select between "Preferences", "About" and "Software update":
 - **Preferences:** Use the drop-down menus to define the preferred unit settings for each parameter. The software needs to be restarted to activate any changes in the settings.
 - **About:** This field gives information about the currently installed software and firmware versions. It allows for starting a firmware update and creating a diagnostics package.
 - **Software Update:** Use this function to download software updates. If connected to the internet, the PC displays an alert '!' in the quick access toolbar when a new software version is released.
- **Quick Access Toolbar (2):** Click this button to create a new experiment.
- **Experiment Menu (4):** The *experiment menu* is already expanded when starting the software. In the *experiment menu*, new experiments can be generated. The *experiment menu* consists of the "Experiment Name", which is typically best defined by the date of experiment and the name of the sample. In the field "User", the user can enter his or her name. The "Location" folder can be changed by clicking on the <...> button. The "Comments" field can be used for any entries concerning the experiment, e.g. information on sample preparation, availability of nitrogen purge or other information. The experiment menu can also be folded and unfolded by clicking on the blue tab "Experiment". A new experiment can be opened by a click on the button <Set up new Experiment> in the *quick access toolbar*. A new experiment must be defined to enable the software to work with the instrument (refer to chapter 5).
- **Current Device Configuration (21):** The *current device configuration* displays the information on all measured parameters in one single

view.

- **Measurement Types:** You can select between six different measurement types:
 - **Dosage (8), not available with SurPASS 3 Eco:** Use the dosage function to either adjust the pH value of the electrolyte solution to a specific pH using one of the titration units, or to dispense a specific volume of liquid by the titration units.
 - **Rinse (9):** Use the rinse function to fill the measuring cell and to rinse the sample with electrolyte solution before starting a zeta potential measurement.
 - **Zeta Potential (10):** Use the zeta potential function to start a zeta potential measurement at the given electrolyte composition.
 - **Isoelectric Point (11), not available with SurPASS 3 Eco:** Use the isoelectric point measurement mode to determine the isoelectric point fully automatically. For each pH step max. 4 zeta cycles are selectable. The start and stop pressure defined in zeta potential are used for the measurement. The start and stop pressure defined in the rinse are used for the rinse cycles.
 - **pH Scan (12), not available with SurPASS 3 Eco:** The pH scan allows to perform a pH dependent zeta potential measurement until a user-defined end pH value is achieved. For each pH step max. 4 zeta cycles are selectable. The start and stop pressure defined in zeta potential are used for the measurement. The start and stop pressure defined in the rinse are used for the rinse cycles.
 - **Adsorption Kinetics (13), not available with SurPASS 3 Eco and only accessible if external pressure supply is activated:** The adsorption kinetics measurement type allows to monitor fast and slow adsorption processes. "Fast adsorption" stops after 3-4 cycles, depending on the gap height. "Slow adsorption" is used for adsorption processes where a longer measurement time is needed.
- **Start Button (20):** Use the <start button> to start the selected measurement type.
- **Measurement View (5):** Select the *measurement view* to monitor the performance of the pressure ramps during measurement or to control the gap adjustment by using the time during rinse cycles.

- **Result View (6):** Select the *result view* to display the result of your zeta potential measurement.
- **Clean (22):** A click on <Clean> opens the *cleaning procedures* with the following functions:
 - **Empty the system:** The task is used for pumping air through the measuring cell and thus removing any liquid after a measurement has finished.



WARNING

Risk of injury

This process must be started with mounted measuring cell with Gap Height/Permeability Index adjusted. If you start the empty process with open gap or the cleaning device mounted, the liquid will splash out of the beaker cover.

- **Clean the system:** This task is used for cleaning the electrolyte system with the cleaning device mounted.
- **Fill the titration units (not available with SurPASS 3 Eco):** This task is used for filling the hoses of the titration unit with acid and base.
- **External Pressure Supply (19):** Click on the symbol to activate or deactivate the external pressure supply and thus the accessibility of the adsorption kinetics feature. The status of the external pressure supply is displayed as green or gray sign.
- **Maintenance (23):** A click on <Maintenance> allows to select between:
 - **Conductivity Calibration:** The conductivity calibration shall be performed once a month. A message will be highlighted when the calibration is due.
 - **pH Calibration:** The pH calibration shall be performed once a week. A message will be highlighted when the calibration is due.
 - **Liquid Declaration (not available with SurPASS 3 Eco):** Assign acid and base to syringe pumps A and B of the titration unit for a correct functioning of measurement types Isoelectric Point and pH Scan.
- **SurPASS 3 Viewer (3):** Click on the <SurPASS 3 Viewer> symbol to open a new SurPASS 3 viewer instance or to bring the viewer to the foreground.

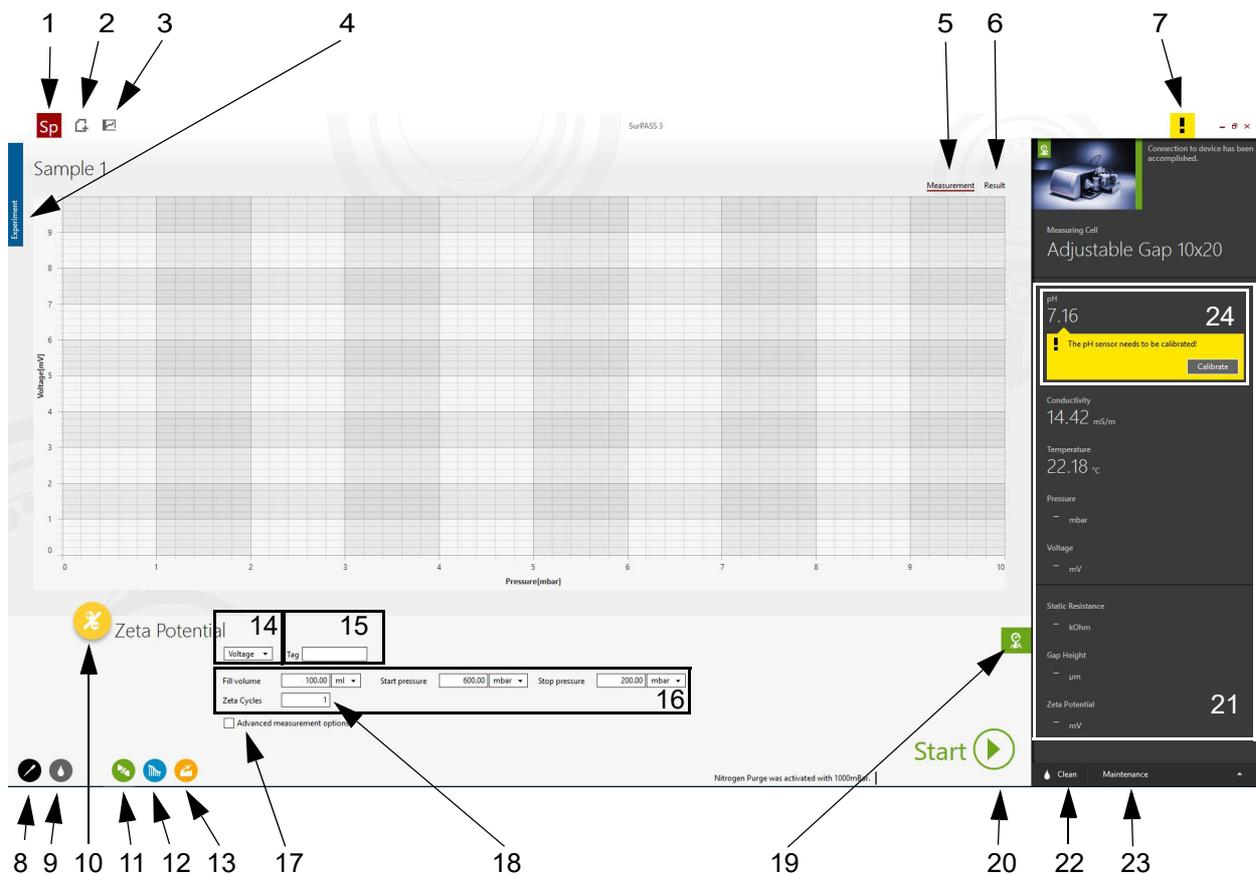


Fig. 19: Graphical user interface of the SurPASS 3 software

- | | |
|---|---|
| <ul style="list-style-type: none"> 1 Application button 2 Quick Access Toolbar for starting a new experiment 3 SurPASS 3 viewer 4 Experiment menu 5 Measurement view 6 Result view 7 Device status 8 Dosage 9 Rinse 10 Zeta Potential 11 Isoelectric Point 12 pH Scan | <ul style="list-style-type: none"> 13 Adsorption Kinetics 14 Selection between voltage and current measurement 15 Tag to identify measurement file 16 Measurement parameters 17 Advanced measurement options 18 Number of measuring cycles 19 Status of the external pressure supply 20 Start button 21 Current device configuration 22 Clean tasks (Empty, Clean and Fill) 23 Maintenance 24 pH calibration reminder |
|---|---|

3.10.3 Export of Data

All measured data are automatically saved in *.xlsx- and *.csv-files. By default result files are saved in the folder `C:\Users\username\Documents\SurPASS3\Author\Experiment` (refer also to chapter 5.6).

TIP: The content of the respective *.xlsx- and *.csv-file is identical, except for the unit settings (customized unit settings in *.xlsx-file, SI units in *.csv-file).

The *.csv-file is needed by the SurPASS 3 viewer and must not be deleted.

SurPASS 3 software serves primarily for the purpose of instrument operation. Measurement files can be re-opened in the SurPASS 3 software, and are then displayed in the SurPASS 3 viewer. A spreadsheet software needs to be used for evaluation of measured data.

4 Putting SurPASS 3 into Operation

4.1 pH Electrode Calibration

The calibration of the pH electrode requires three buffer standard solutions of known pH value. The buffer standard solutions for calibrating the pH electrode should equally cover the complete measuring range (e.g. pH 3, pH 7, pH 10). To ensure an accurate calibration and pH measurement it is important to use fresh buffer solutions for the calibration.

NOTICE

- It is recommended to calibrate the pH electrode once a week. A message will be highlighted when the calibration is due.
- If not in use store the pH electrode in 3 mol/l KCl solution and place it in the storage container of the accessory kit for liquid handling (pos. 5 in Fig. 17).

1. Click on <Maintenance> and open the task "pH calibration".

TIP: If the message for the need for pH calibration is displayed, a click on the message guides you to the task "pH calibration".

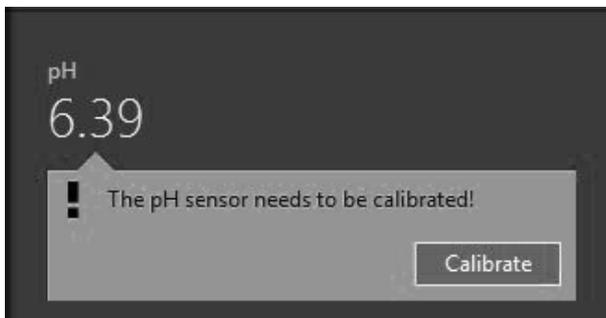


Fig. 20: pH calibration reminder

- Click on <Begin calibration>.
- Clean the pH electrode with ultrapure water and insert it into the first buffer solution.
- Enter the pH value of this first buffer solution in the first number field. The measured voltage and pH value will be displayed. Click on the small triangle next to "Start" to start the pH cali-

bration.

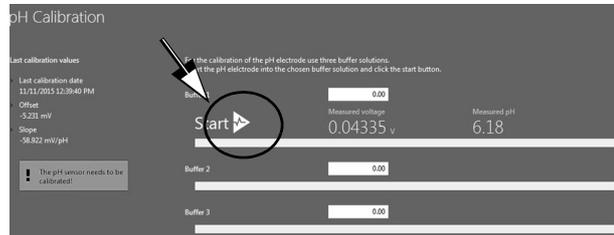


Fig. 21: pH calibration

- The progress bar indicates the progress of the measurement. Wait until it has finished, then clean the pH electrode with water and place it into the next buffer solution. Enter its pH value into the second number field and click on the <Start> button to continue the pH calibration.
- Proceed accordingly with the third buffer solution.

NOTICE

At 25 °C the following values for the pH calibration parameters Slope and Offset are acceptable:
Slope: -59.16 mV/pH ± 4 % (-56.8 mV/pH to -61.5 mV/pH)
Offset: ± 30 mV

- If the pH calibration parameters are accepted by the software, click on <Accept> to save them. A green checkmark will indicate a successful calibration.
- If any of the values is out of range, click on <New calibration> to repeat the pH calibration, or click on <Close> to repeat the pH calibration at a later time.

Possible reasons for an unsuccessful pH calibration are:

- an invalid pH of one of the buffer standard solutions (consider a temperature effect on pH),
- a drift of the pH electrode or
- a damage of the pH electrode connector or cable.

4.2 Conductivity Electrode Calibration

The calibration of the conductivity electrode is a single point calibration and requires a 0.1 mol/l KCl solution or a conductivity standard solution.

The conductivity electrode delivered with the SurPASS 3 contains a temperature sensor. The temperature measurement by this sensor is used for both conductivity calibration and zeta potential evaluation.

NOTICE

- It is recommended to calibrate the conductivity electrode once a month. A message will be highlighted when the calibration is due.
- If not in use store the conductivity electrode in dry conditions after cleaning and place it in the storage container of the accessory kit for liquid handling (pos. 4 in Fig. 17).

1. Click on <Maintenance> and open the task "Conductivity calibration".

TIP: If the message for the need for conductivity calibration is displayed, a click on the message guides you to the task "conductivity calibration".

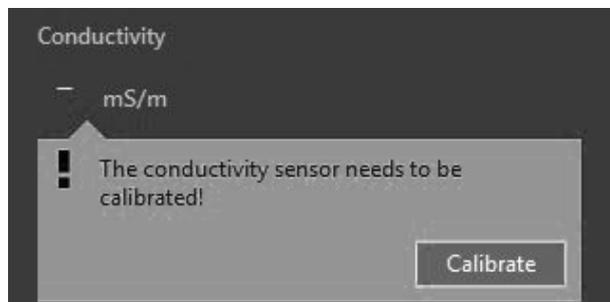


Fig. 22: Conductivity calibration reminder

2. Click on <Begin calibration>.
3. Choose a 0.1 mol/l solution of KCl or a conductivity standard for conductivity calibration.
4. Prepare 200 ml of the 0.1 mol/l KCl solution or conductivity standard solution in the 250 ml glass beaker.

4.2.1 Conductivity Calibration using 0.1 mol/l KCl

1. Select "0.1 mol/l KCl" as calibration liquid.

2. Clean the conductivity electrode with ultrapure water and insert it into the 250 ml glass beaker filled with 0.1 mol/l KCl solution until the hole in the shaft is immersed completely. The measured conductivity and temperature will be displayed.
3. Click on the small triangle next to "Start" to start the conductivity calibration.

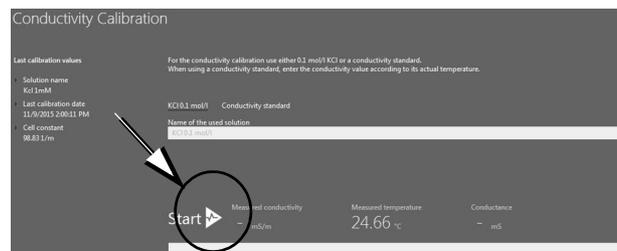


Fig. 23: Conductivity calibration

4. The progress bar indicates the progress of the measurement. Wait until it has finished.
5. If the cell constant is accepted by the software, click on <Accept> to save it. A green checkmark will indicate a successful calibration.
6. If the cell constant is unacceptable, click on the small triangle to start the calibration again, or click on <Close> to repeat the conductivity calibration at a later time.

TIP: The conductivity electrode supplied with the SurPASS 3 instrument has a cell constant of $47.5 \text{ m}^{-1} \pm 5.25 \% (45 \text{ m}^{-1} \dots 50 \text{ m}^{-1})$. A persistent deviation by $> 5.25 \%$ indicates the needs for exchanging the conductivity electrode.

4.2.2 Conductivity Calibration using a Conductivity Standard

1. When using a conductivity standard, select "conductivity standard" and enter the name of the conductivity standard and its conductivity at the given temperature as stated on the bottle.
2. Proceed with steps 2-6 in chapter 4.2.1.

TIP: The conductivity electrode supplied with the SurPASS 3 instrument has a cell constant of $47.5 \text{ m}^{-1} \pm 5.25 \% (45 \text{ m}^{-1} \dots 50 \text{ m}^{-1})$. A persistent deviation by $> 5.25 \%$ indicates the needs for exchanging the conductivity electrode.

TIP: Watch the correct unit for conductivity when using a conductivity standard.

4.3 Filling of the Titration Unit (not applicable to SurPASS 3 Eco)

The SurPASS 3 instrument with titration unit contains two syringes with a volume of 1.0 ml for dispensing acid and base to change the pH of the electrolyte solution.

The hoses of the titration unit have to be filled before starting a pH titration. If different titration liquids are in use, the hoses have to be rinsed after exchanging the titration liquid. Make sure the titration hoses are mounted properly before proceeding with filling the titration units.

1. In order to assign the liquids in use to the respective syringe pump, select <Maintenance> and "Liquid Declaration".
2. Select "Acid", "Base" or "Other" for each syringe pump and save these settings.

TIP: "Acid" or "Base" needs to be selected in order to enable isoelectric point or pH scans, respectively.

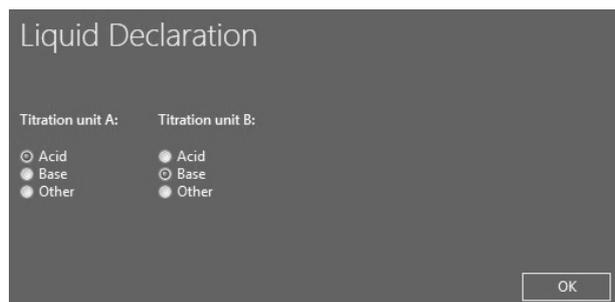


Fig. 24: Liquid declaration

3. In order to fill the titration unit, select <Clean> and "Fill the titration units".
4. Click on <Start> to start the filling of the titration units. The system uses a volume of 4 ml for each syringe pump to ensure proper filling of the hoses.

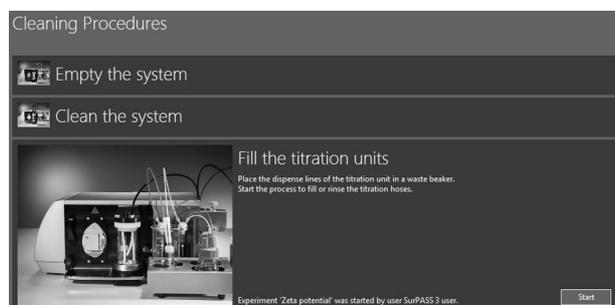


Fig. 25: Fill the titration units

TIP: The same procedure can be used for cleaning and emptying the titration unit. Use ultrapure water for cleaning and put the dispense line in a waste beaker. Remove the uptake line from the liquid reservoir to empty the titration unit by performing fill cycles with air.



WARNING

Risk of injury

A malfunction of the titration unit may introduce a high concentration of hazardous liquids into the SurPASS 3 electrolyte circuit. Wear protective clothing during cleaning.



CAUTION

Make sure to have the titration burette tips removed from the 250 ml beaker when switching on the SurPASS 3 instrument. During initialization the syringes move and dispense of titration liquid into the electrolyte beaker may take place unintentionally.

4.4 SurPASS 3 Verification Measurement with Cotton Fabrics

Follow these instructions in order to perform a verification measurement using a standard cotton fabric (ISO 105-F09) on your SurPASS 3 instrument. Meeting the specified range of the zeta potential of this cotton fabric confirms that the SurPASS 3 is in a reliable working condition.

4.4.1 Chemicals and Accessories

In order to perform the verification measurement, you need:

1. the accessory for SurPASS 3 verification measurement (cat.no. 181830) incl. a set of standard fabrics for verification measurement (cat.no. 182927)



Fig. 26: Scope of supply of the accessory for SurPASS 3 verification measurement

2. KCl, p.A.
3. ultrapure water (ASTM I grade)
4. precision balance

4.4.2 Preparation of the 0.01 mol/l KCl Standard Solution for the Verification Measurement with Ag/AgCl Electrodes

1. Rinse the 250 ml volumetric flask that is contained in the accessory for SurPASS 3 verification measurement with ultrapure water.
2. Add the amount of KCl required to prepare the 0.01 mol/l standard solution, *i.e.* 0.186 g for 250 ml.

Volume of the volumetric flask [ml]	Amount of KCl [g]	Amount of KCl [mg]
250 ^a	0.186	186
500	0.373	373
1000	0.746	746

^a Contained in the accessory for SurPASS 3 verification measurement

3. Transfer the KCl salt into the volumetric flask. Make sure that the complete amount of KCl is transferred.
4. Fill the volumetric flask with ultrapure water exactly up to the ring marker.

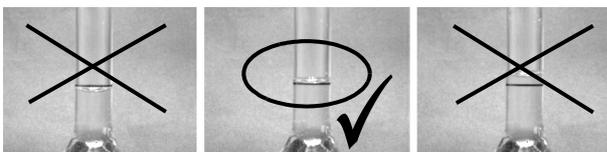


Fig. 27: Preparation of the measuring liquid

5. Close the volumetric flask and tilt it several times to guarantee a homogeneous mixing of the salt with the ultrapure water.

4.4.3 Preparation of the 0.001 mol/l KCl Standard Solution for the Verification Measurement with Gold Electrodes

1. Rinse the 250 ml volumetric flask that is contained in the accessory for SurPASS 3 verification measurement with ultrapure water.
2. Add the amount of KCl required to prepare the 0.001 mol/l standard solution.

Volume of the volumetric flask [ml]	Amount of KCl [g]	Amount of KCl [mg]
250 ^a	0.0186	18.6
500	0.0373	37.3
1000	0.0746	74.6

^a Contained in the accessory for SurPASS 3 verification measurement

3. Transfer the KCl salt into the volumetric flask. Make sure that the complete amount of KCl is transferred.
4. Fill the volumetric flask with ultrapure water exactly up to the ring marker.

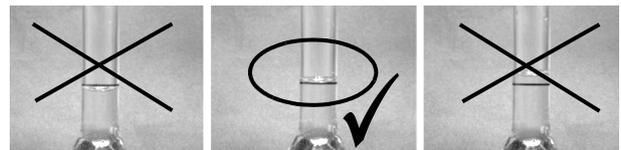


Fig. 28: Preparation of the measuring liquid

5. Close the volumetric flask and tilt it several times to guarantee a homogenous mixing of the salt with the ultrapure water.

4.4.4 Mounting the Standard Cotton Fabric

1. Take one standard cotton fabric and fold it four times. You get a square sample of 1 cm x 1 cm with 16 cotton fabric layers.

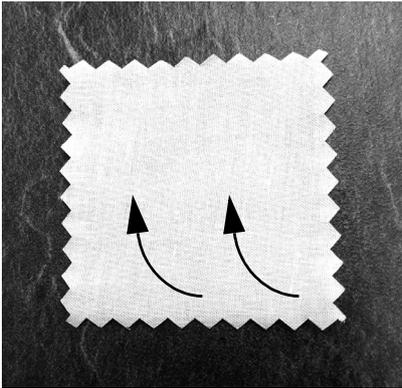


Fig. 29: Folding step 1

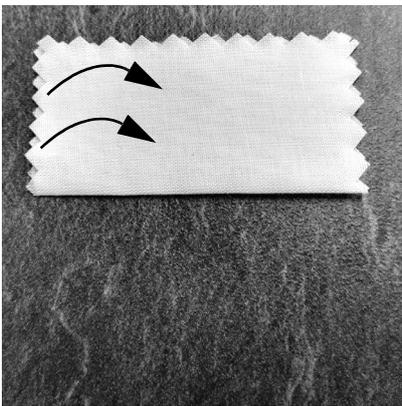


Fig. 30: Folding step 2

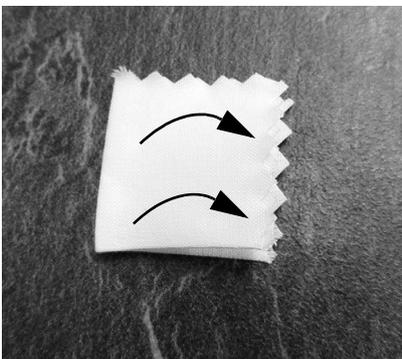


Fig. 31: Folding step 3

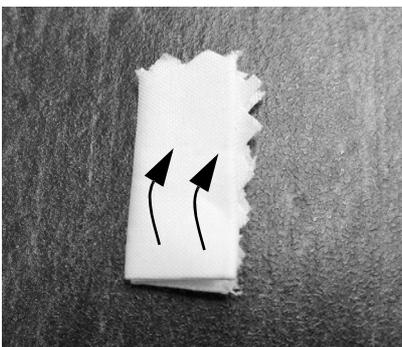


Fig. 32: Folding step 4



Fig. 33: Final size (1 cm x 1 cm) of the folded cotton fabric

2. Mount the folded cotton fabric into the clean sample holder of the SurPASS 3 verification device. Position the fabric in the center of the sample holder.

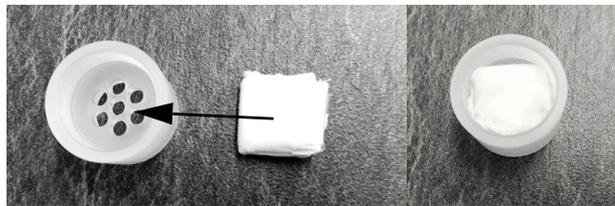


Fig. 34: Mounting the cotton fabric

3. Make sure that no fibers are blocking the thread when closing the sample holder.
4. Use the two mounting tools to tighten the sample holder completely, so that no gap is visible between the two halves.

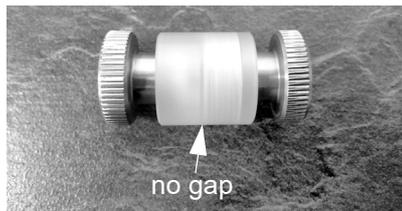


Fig. 35: Closed sample holder with mounting tools

5. Mount the sample holder into the SurPASS 3 verification device. Fix the screws by using the Torx T20 screwdriver with torque control.
6. Mount the SurPASS 3 verification device on the instrument as described in chapter 5.1.
7. Transfer 200 ml of the 0.01 mol/l KCl (or 0.001 mol/l KCl) volumetric electrolyte solution into the 250 ml glass beaker.
8. The conductivity of the 0.01 mol/l KCl (or 0.001 mol/l KCl) standard solution has to match the tabulated data. A deviation of < 5 % is acceptable.

Table 4-1: Conductivity data

temperature in °C	0.01 mol/l KCl solution	
	conductivity in mS/m	conductivity in $\mu\text{S}/\text{cm}$
20	127.6	1276
21	130.4	1304
22	133.1	1331
23	135.8	1358
24	138.6	1386
25	141.3	1413
26	144.1	1441
27	146.8	1468
28	149.6	1496
29	152.4	1524
30	155.0	1550

Source: Handbook of Chemistry and Physics

9. The pH of the 0.01 mol/l KCl standard solution has to fit within the limits of pH 5.5 - pH 6.5.

4.4.5 Verification Measurement with Standard Cotton Fabric

1. Start a new experiment in the SurPASS 3 software as described in chapter 5.2.

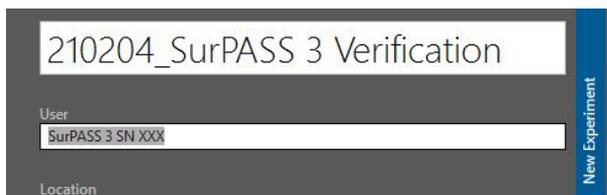


Fig. 36: Experiment for SurPASS 3 verification measurement

2. Start 5 rinse cycles with the default parameters:
 - Fill volume: 100 ml
 - Start pressure: 600 mbar
 - Stop pressure: 200 mbar
 - Rinse cycles: 5



Fig. 37: Rinse steps

3. Verify that the permeability index is in the range of 100 ... 130.
4. Start 5 Zeta Potential measurements with the default parameters as described in chapter 5.3.1:
 - Fill volume: 100 ml
 - Start pressure: 600 mbar
 - Stop pressure: 200 mbar
 - Zeta cycles: 5
 - Activate "Advanced measurement options"
 - Upper pressure limit: 450 mbar

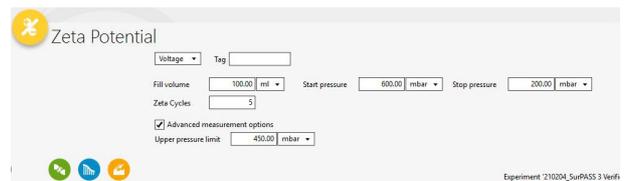


Fig. 38: Zeta potential measurements

4.4.6 Verification Results for Ag/AgCl Electrodes

Use the results of measurements no. 3, no. 4 and no. 5 to evaluate the zeta potential of the standard cotton fabric. The expected value for the zeta potential is -19 mV. A deviation of $\pm 10\%$ is acceptable.

The corresponding pressure ramp data (streaming potential vs. pressure difference) shall be linear and overlapping.



Fig. 39: Pressure ramps

TIP: Refer to the reference guide D85IB001EN for more information, tips and tricks in case the verification measurement did not yield the expected result.

4.4.7 Verification Results for Gold Electrodes

Use the results of measurements no. 3, no. 4 and no. 5 to evaluate the zeta potential of the standard cotton fabric. The expected value for the zeta potential is $-15 \text{ mV} \pm 2 \text{ mV}$.

4 Putting SurPASS 3 into Operation

The corresponding pressure ramp data (streaming potential vs. pressure difference) should be linear and overlapping.



Fig. 40: Pressure ramps with gold electrodes

5 Performing a Measurement

NOTICE

The SurPASS 3 was tested and cleaned before shipment. It is not necessary to rinse or clean the instrument before starting the first measurement after installation.

NOTICE

Make sure to use ultrapure water (ASTM I) and high purity chemicals as described in chapter 2.4 before starting the first measurement.

In order to perform a successful zeta potential analysis with SurPASS 3, the following procedures have to be performed in the given order:

1. Mount the sample in a suitable measuring cell (refer to the reference guide D85IB001EN).
2. Mount the measuring cell on SurPASS 3 (see chapter 5.1).
3. Define a new experiment in the software and rinse the system with electrolyte solution (see chapter 5.2).
4. Adjust the gap height if one of the Adjustable Gap Cells or the Clamping Cell is used (see chapter 5.2.1).
Adjust the permeability index if the Cylindrical Cell is used (see chapter 5.2.2).
5. Define the parameters for either a zeta potential analysis, isoelectric point determination or pH scan and start the measurement (see chapter 5.3).
6. After finishing the measurement, empty the system (see chapter 5.5.1).
7. Remove the measuring cell with the sample and mount the cleaning device (see chapter 5.5.2).
8. Clean the system by rinsing the electrolyte circuit with ultrapure water or an appropriate cleaning solution (see chapter 5.5.3).

5.1 Mounting the Measuring Cells on SurPASS 3

1. Observe the correct positioning of the measuring cell with the arrow always pointing upwards.



Fig. 41: Arrow on the measuring cell pointing upwards

2. Make sure that all parts are dry and that there is no liquid between the electrodes and the measuring cell.
3. Push the measuring cell into the interlock and turn the locking lever on the SurPASS 3 upwards.



Fig. 42: Locking lever



CAUTION

Risk of injury

Watch your fingers when fixing the measuring cell.

4. The SurPASS 3 software will recognize that a measuring cell is connected and the name of the measuring cell will be displayed in the *current device configuration*.

5.2 Rinsing the System

NOTICE

- Make sure the complete electrolyte circuit and the measuring cell were rinsed with ultra-pure water (ASTM I) before starting a measurement.
- Make sure the pH and conductivity electrodes are calibrated.

1. Check the proper connection of the external electrolyte hoses (see chapter 3.5).

NOTICE

The water used for rinsing and for preparing the electrolyte solution for zeta potential measurement must conform with the American Society for Testing and Materials standard ASTM I.

NOTICE

Do not clean the glass bottles used with SurPASS 3 in the dishwasher. Residues of detergents are difficult to remove from the SurPASS 3 electrolyte circuit and can falsify the zeta potential results.

2. Put the magnetic stir bar and fill 200 ml of electrolyte solution into the glass beaker.
3. If required start purging the electrolyte solution with nitrogen. For details on the requirements for nitrogen supply refer to appendix C. Make sure that the external pressure is activated in the software as indicated by a green external pressure supply sign (refer to chapter 3.10.2).
4. Create a new experiment in the SurPASS 3 software by filling out the fields "Experiment Name" and "User". The file "location" can be altered by clicking on <...> and selecting a different folder. Any comments on the experiment can be entered and the new experiment is started by clicking on <Set up new Experiment>.

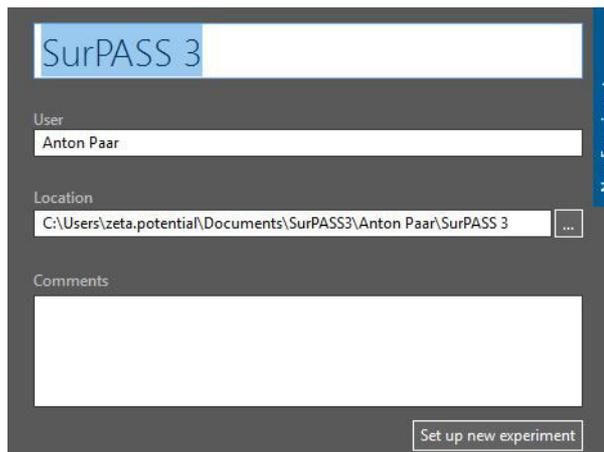


Fig. 43: New experiment

TIP: Typically, the name of the sample is an appropriate name for the experiment.

TIP: The new experiment menu is already open when starting the software. If you need to open the menu manually, click on the <Ste up new Experiment> button in the quick access toolbar.

5. Select the task "Rinse". Use the default values of 100 ml as fill volume, 600 mbar as start pressure, 200 mbar as stop pressure and 3 cycles.



Fig. 44: Rinse cycle

TIP: Make sure that nitrogen is available and that the external pressure is activated in the software (refer to chapter 3.10.2) if the start pressure needs to be set higher than 1500 mbar.

6. Click on <Start>.
7. Double-click into the graph for full display of the data. Follow the performance of the pressure ramps ramp (pressure vs. time) in the *measurement view*.
8. If necessary adjust the gap height or the permeability index during the rinse cycles.

TIP: If the streaming resistance (refer to chapter 5.2.1) is extremely high, residues of air are remaining in the measuring cell. To get rid of the air, open the gap by using the adjustment knob (for Adjustable Gap Cell and Clamping Cell) during the rinse cycle and then close it again.

TIP: The defined values for start and stop pressure are also used for the rinsing cycles during the "Isoelectric Point" and "pH Scan" measurement

types.

5.2.1 Adjusting the Gap between Planar Sample Surfaces

1. Monitor the pressure drop over time which is displayed continuously in the *current device configuration*. Adjust the gap height by rotating the adjustment knob.

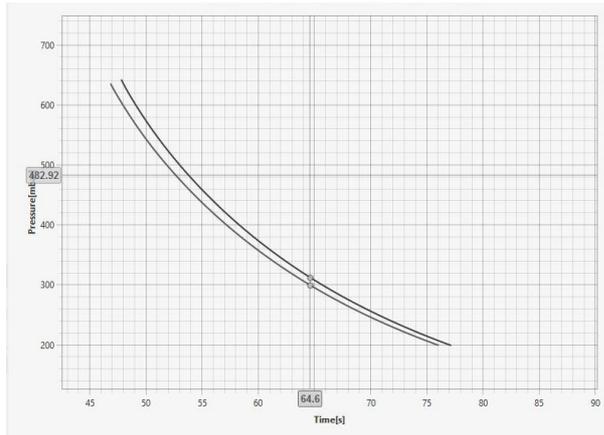


Fig. 45: Rinse cycles

TIP: With the standard measurement parameters, the end pressure of 200 mbar should be reached in approx. 65 to 75 seconds so that your alignment is in the range of 80 to 120 μm .

2. Monitor the gap height which is displayed in the *current device configuration* after each rinse cycle has finished.

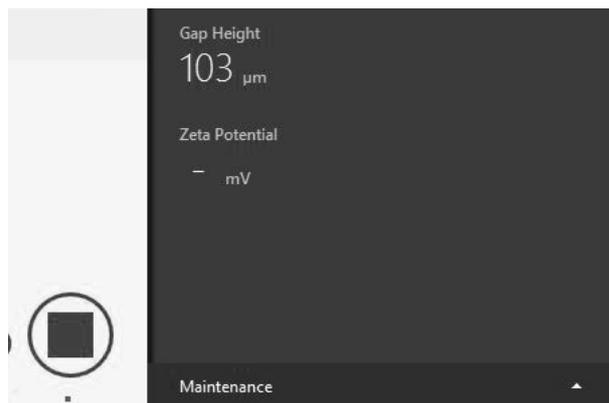


Fig. 46: Gap height

3. According to this gap height,
 - a. use the adjustment knob to close/open the gap carefully to achieve the target gap of 100 μm (1 unit of the scale on the adjustment knob equals 25 μm). If necessary start additional rinse cycles.

- b. close the adjustment knob by min. 300 μm , if no gap is displayed
- c. open the adjustment knob if the pressure displayed in the *current device configuration* does hardly decrease (streaming resistance is extremely high).

4. Repeat steps 2-3 until the gap height is adjusted to 100 μm .

5.2.2 Adjusting the Permeability Index of Sample Plugs in the Cylindrical Cell

1. Monitor the permeability index which is displayed in the *current device configuration* after the rinse cycle has finished.
2. Adjust the permeability index approximately to 100 by turning the micrometer screw.

5.3 Starting a Zeta Potential Measurement

The routines "Rinse" (chapter 5.2), "Zeta Potential" (chapter 5.3.1), "Isoelectric Point" (chapter 5.3.2) and "pH Scan" (chapter 5.3.3) record pressure ramps. Fill volume, start pressure and stop pressure can be defined for each of these routines. The default settings (Fill volume 100 ml, Start pressure...) can be used for most analyses.

5.3.1 Zeta Potential

1. Select the measurement type "Zeta Potential" to perform a single or a series of zeta potential measurements at defined electrolyte composition.

TIP: If you need to change the pH value of your electrolyte solution, first select the task "Dosage" to either add acid or base until a certain pH value is reached or to add a certain volume of acid or base. Make sure to insert the corresponding burette tip into the electrolyte beaker before clicking on <Start>. Perform at least two rinse cycles (see chapter 5.2) before starting a zeta potential measurement after changing the pH value.

2. Choose one (1) cycle if you wish to perform a single zeta potential measurement, or type in a higher number of measurements (up to 99) for a series of zeta potential measurements at same electrolyte composition.
3. Click on the <Start> button.

5 Performing a Measurement



Fig. 47: Zeta potential measurement

4. Monitor the performance of the pressure ramps in the *measurement view* during measurement by a double-click in the graph.

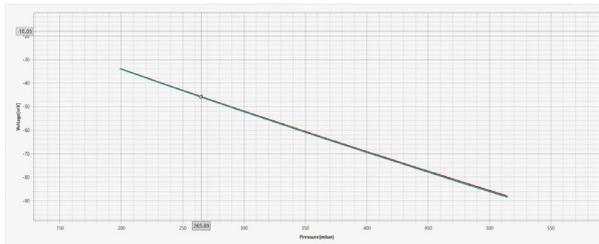


Fig. 48: Pressure ramps (measurement view)

5. The performance of the pressure ramps is very important, since the software is doing a linear fit through the recorded streaming current / streaming potential values.
If necessary use the Advanced measurement options to define the upper pressure limit.



Fig. 49: Advanced measurement options

Check the box "Advanced measurement options" and type in the upper pressure where the recorded pressure ramp still appears linear, e.g. 450 mbar.

6. Find the zeta potential result in the *current device configuration* and in the *result view* after the measurement has finished.

5.3.2 Isoelectric Point (not applicable to SurPASS 3 Eco)



WARNING

Hazardous Liquids

The volume of electrolyte solution in the beaker must not exceed the 200 ml mark of the beaker when starting an Isoelectric Point or pH Scan to avoid an overflow of liquid during the course of the analysis.

1. Select the measurement type "Isoelectric Point" for the most efficient way to determine the isoelectric point of your sample.
2. Type in a number of rinse cycles which the instrument shall execute after the pH value was adjusted and before performing another zeta potential measurement at changed electrolyte composition. The start and stop pressure, defined in "Rinse", are used for these rinse cycles (refer to chapter 5.2).

TIP: Three rinse cycles are sufficient for most samples. If you observe from the measurement view that the pressure ramps are not overlapping after three cycles, stop the measurement and use a higher number of cycles to ensure equilibrium conditions for correct zeta potential determination.

3. Type in a number of zeta cycles which the instrument shall execute at the adjusted pH value. A max. of 4 repetitions are possible.
4. For the measurement the values for start and stop pressure defined in "Zeta Potential" are used.
5. The pH increment is set by default to 0.3 pH units and is adjustable up to 2 pH units.
6. Wipe off the burette tip of both titration units and place them into the electrolyte beaker. Make sure the titration hoses are filled with the respective titration liquids (refer to chapter 4.3).
7. Click on the <Start> button.



Fig. 50: Isoelectric point determination

8. Monitor the performance of the pressure ramps in the *measurement view* during measurement by a double-click in the graph.
9. Find the zeta potential vs. pH dependence and the isoelectric point in the *result view* after the measurement has finished.

5.3.3 pH Scan (not applicable to SurPASS 3 Eco)



WARNING

Hazardous Liquids

The volume of electrolyte solution in the beaker must not exceed the 200 ml mark of the beaker when starting an Isoelectric Point or pH Scan to avoid an overflow of liquid during the course of the analysis.

1. Select the measurement type "pH scan" for zeta potential determination in a defined pH range.
2. Type in a number of rinse cycles which the instrument shall execute after the pH value was adjusted and before performing another zeta potential measurement at changed electrolyte composition. The start and stop pressure, defined in "Rinse", are used for these rinse cycles (refer to chapter 5.2).

TIP: *Three rinse cycles are sufficient for most samples. If you observe from the measurement view that the pressure ramps are not overlapping after three cycles, you may want to stop the measurement and use a higher number of cycles to ensure equilibrium conditions for correct zeta potential determination.*

3. Type in a number of zeta cycles which the instrument shall execute at the adjusted pH value. A max. of 4 repetitions are possible.
4. For the measurement the values for start and stop pressure defined in "Zeta Potential" are used.
5. Type in the end pH value where you want the measurement to stop. The instrument will perform zeta potential measurements starting from the pH value of the electrolyte solution until the end pH value was passed. The pH increment is set by default to 0.3 pH units and is adjustable up to 2 pH units.
6. Wipe off the burette tip of the titration unit which is filled with acid or base, depending on the just defined end pH value, and place it into the electrolyte beaker. Make sure the titration hoses are filled with the respective titration liquid (refer to chapter 4.3).

7. Click on the <Start> button.



Fig. 51: pH scan

8. Monitor the performance of the pressure ramps in the *measurement view* during measurement by a double-click in the graph.
9. Find the zeta potential vs. pH dependence in the *result view* during and after the measurement has finished.

5.4 Adsorption Kinetics Measurement (not applicable to SurPASS 3 Eco)

Please refer to the reference guide D85IB001EN for more information about adsorption kinetics measurement.

5.5 End of Measurement

A measurement has finished when the <Stop> button turns into a <Start> button again.

5.5.1 Empty the System

1. Click on <Clean> and select the task "Empty the system" in the *cleaning procedures*.



Fig. 52: Empty the system

2. Leave both the outlet and the inlet hoses in the beaker with the electrolyte solution.

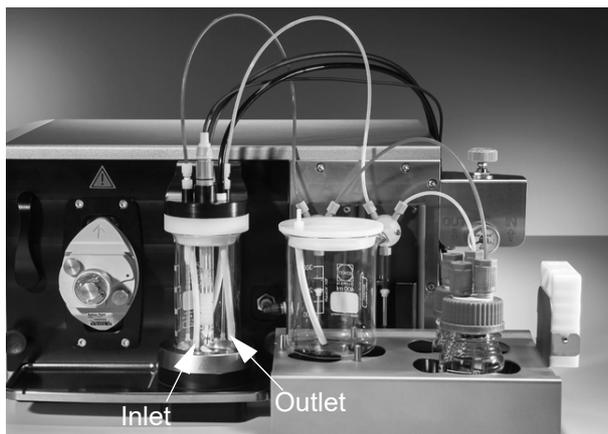


Fig. 53: System configuration while emptying the system

3. Click on <Start> to pump air through the measuring cell and thus remove all liquid from the electrolyte circuit.

NOTICE

Avoid starting an empty process while the cleaning device or a cell with too large gap height is mounted as the measurement solution will splash out of the beaker.



Fig. 54: Turn the locking lever

2. Pull out the measuring cell for sample removal.



Fig. 55: Take off the measuring cell

5.5.2 Removing the Measuring Cell



WARNING

Risk of injury

Do not remove the measuring cell while the instrument is in operation.



CAUTION

Hazardous liquids

Always empty the electrolyte circuit of SurPASS 3 before removing the measuring cell.

1. Turn the locking lever on the front panel to the left.

5.5.3 Clean the System

1. Lift the cover or the 250 ml electrolyte beaker and connect the outlet hose (on the right side) to the beaker cover of the waste beaker.
2. Fill 200 ml of ultrapure water or cleaning solution into the 250 ml beaker.

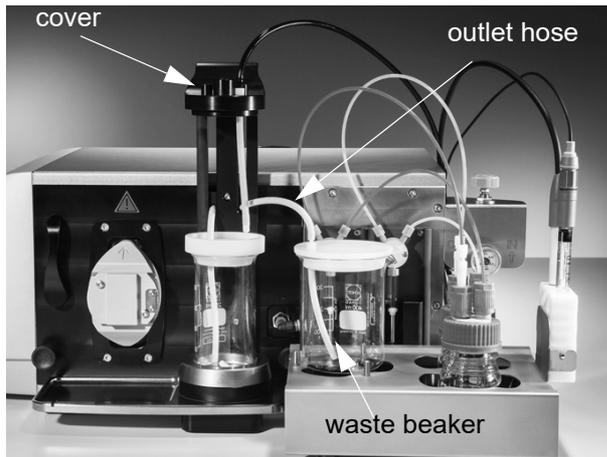


Fig. 56: System configuration while cleaning the system



WARNING

Risk of injury

Do not use inflammable, toxic or explosive liquids for SurPASS 3 operation (see chapter 1.1).

3. Click on <Clean> and select the task "Clean the system" in the *cleaning procedures*.

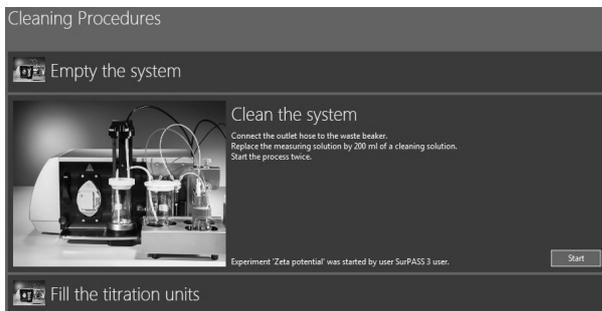


Fig. 57: Cleaning the system

4. Mount either the adapter for cleaning or the measuring cell on the SurPASS 3 instrument (see chapter 5.1).
5. Start the clean cycle by clicking on <Start>.
6. If necessary repeat steps 1-5 with a fresh cleaning solution.

5.5.4 Switching off the Instrument

NOTICE

Before switching off the SurPASS 3 instrument, make sure the electrolyte system is clean and empty (refer to chapter 5.5.3).

1. Switch off the nitrogen supply in your laboratory.
2. Switch off the instrument by using the power switch on the rear side of the SurPASS 3 instrument.
3. Store the pH electrode in 3 mol/l KCl solution and place it in the storage position of the liquid handling kit (Pos. 5 in Fig. 17).
4. Place the conductivity electrode in dry condition into its storage position of the liquid handling kit (Pos. 4 in Fig. 17).



WARNING

Risk of injury

When switching off the SurPASS 3 software and instrument independently of the nitrogen supply, the hose connecting the SurPASS 3 instrument and your in-house gas supply (hose (5) in Fig. 18) will remain pressurized.

It needs to be depressurized in case you intend to disconnect it (see chapter 6.9).

TIP: If you first switch off the nitrogen supply but keep the external gas supply activated in the software and the instrument switched on, a strange noise will indicate that the system is depressurized. Switch off the instrument to stop the noise.

5.6 Data Analysis

5.6.1 Export of Measured Data

After a measurement has finished, all data is saved in the folder which was defined as file location upon creating the experiment. By default, this is:

C:\Users\username\Documents\SurPASS3\Author\Experiment name (for Windows 7)

C:\My computer\Documents\SurPASS3\Author\Experiment name (for Windows 8)

In this folder, you will find:

- for an individual zeta potential measurement:
 - *.xlsx- and *.csv-files with the results for each

5 Performing a Measurement

zeta potential measurement

- a subfolder containing *.xlsx- and *.csv-files of the raw data for each measurement
 - the*.surf-file for display of the measured data in the SurPASS 3 viewer (see reference guide D85IB001EN).
- for an isoelectric point determination:
 - *.xlsx- and *.csv-files of the complete IEP scan
 - a subfolder containing *.xlsx- and *.csv-files of the raw data for each zeta potential measurement
 - the*.surf-file for display of the measured data in the SurPASS 3 viewer (see reference guide D85IB001EN).
- for a pH scan:
 - *.xlsx- and *.csv-files of the complete pH scan
 - a subfolder containing *.xlsx- and *.csv-files of the raw data for each zeta potential measurement
 - the*.surf-file for display of the measured data in the SurPASS 3 viewer (see reference guide D85IB001EN).

TIP: *The content of the respective *.xlsx- and *.csv-file is identical, except for the unit settings (customized unit settings in *.xlsx-file, SI units in *.csv-file). The *.csv-file is needed by the SurPASS 3 viewer and must not be deleted.*

6 Upkeep and Cleaning

This chapter contains information on how to handle the instrument to ensure its smooth long-term operation.

- After a standby > 2 weeks, or after a series of different application measurements, it is recommended to perform the instrument verification as described in section 4.4 SurPASS 3 Verification Measurement with Cotton Fabrics.
- Please take note of section 4.1 and section 4.2 concerning the pH and conductivity electrode calibration:
 - The pH electrode should be calibrated once a week.
 - The conductivity electrode should be calibrated once a month.
- To assure a constant and high accuracy of your measurement, operate and store the instrument under the specified conditions and perform the recommended tasks regularly.
- Ensure that all environmental requirements described in section 1.1 General Safety Instructions (sub-section "Installation and use") as well as in section 2.3 System Requirements and section 2.4 Water Quality and Measurement Conditions are always fulfilled.
- Always handle the instrument with care. Even minor damage may cause serious measurement errors.
- Employ a regular optical check if there is no contamination on the outside of the instrument and perform an effective cleaning routine.



WARNING

Risk of injury

Repair and service activities exceeding the procedures described in this instruction manual may only be carried out by authorized personnel. In case of a malfunction of the SurPASS 3 instrument contact your local Anton Paar representative.

6.1 Software Administration

- Keep the delivered USB stick with the software delivered with the instrument in a safe place.
- Perform an update if the software indicates a

new firmware or software version available in the cloud.

- Make sure that the used computer meets the needed specifications as described in appendix A Technical Data and in section 2.3 System Requirements.

6.2 Upkeep and Cleaning after Standard Operation

- Wipe the instruments cover to keep it free of dust and to avoid chemical damage on the instrument cover surface.
- Perform a clean cycle with ultra-pure water after the measurement operation and before switching off the instrument.
- It is not allowed to operate or clean the instrument with flammable liquids.
- Rinse the pH electrode with ultra-pure water and then put it in the appropriate storage beaker filled with KCL 3 mol/L if the instrument is not in use.
- Rinse the conductivity electrode with ultra-pure water and then put it in the storage beaker of the instrument in dry condition if the instrument is not in use.
- Turn off the nitrogen purge in the software and also the in-house nitrogen supply if not in use.
- Empty the electrolyte and the waste beakers if the instrument is not in use.
- Put the dispense lines for acid and base into the storage position if they are not in use.

6.3 Upkeep and Cleaning after Adsorption Operation

Clean the instrument using ultra-pure water or an appropriate cleaning solution.

To keep the measurement performance an extensive cleaning after the adsorption tests is recommended. Refer to section 6.7 Manual Cleaning of the Electrolyte Circuit.

Note: *It is not allowed to clean the instrument with flammable liquids.*

6.4 Cleaning and Drying the Measuring Cells

- The gaskets and O-rings of all measuring cells need to be checked after a measurement to guarantee trouble-free operation.
- If the cells are not in use, store them disassembled in dry condition.
- Make sure that they were cleaned with ultra-pure water and that all residues of samples are removed.

6.5 Wetted Parts

During the rinse and measurement processes the electrolyte gets in contact with the following materials.

- PharMed BPT tube (fluorinated polymer)
- Silicone tube
- PVDF poly(vinylidene fluoride)
- Borosilicate glass
- SiC Silicon carbide
- FPM/FKM (fluorinated rubber)
- EPDM ethylene propylene diene monomer
- POM-C (Polyoxymethylene Copolymer)

SurPASS 3 adapter for cleaning (cleaning device)

- PP polypropylene

Cylindrical Cell

- PA poly(amide)
- PCTFE poly chloro trifluoro ethylene
- PEEK polyether ether ketone

Adjustable Gap Cell D14/D15

- PPS polyphenylene sulfide

Adjustable Gap Cell 20 x 10 mm

- PEEK GF30 polyether ether ketone

Titration Unit

- ETFE ethylene tetrafluoro ethylene
- PTFE poly(tetrafluoro ethylene)

Clamping Cell

- PEEK GF30 polyether ether ketone

Electrodes

- Ag/AgCl electrodes

NOTICE

Risk of damage.

Do not use the Ag/AgCl electrodes with:

- Ammonia and ammonia generating solutions
- Solutions containing salts of bromide or iodide
- Solutions containing metal ions such as copper.

- alternatively gold electrodes



CAUTION

Risk of damage.

The SurPASS 3 instrument is intended for use with dilute aqueous solutions of inorganic salts, acids and bases.

Before using any additive or a solvent different from water, you have to make sure that all of the materials listed above are chemically compatible.



WARNING

Risk of injury

Do not use inflammable, toxic or explosive liquids for SurPASS 3 operation (see chapter 1.1).

6.6 Exchange of the Electrodes

With SurPASS 3 you may use Ag/AgCl or gold electrodes.

Please refer to the reference guide D85IB001EN for more information about exchanging of the electrodes.

6.7 Manual Cleaning of the Electrolyte Circuit

This chapter contains information on how to clean the electrolyte circuit manually in case of a severe

contamination.

Please refer to the reference guide D85IB001EN for more information about manual cleaning of the electrolyte circuit.

6.8 Removing and Mounting of the Syringes of the Titration Unit (not applicable to SurPASS 3 Eco)

Follow the instructions given in the reference guide D85IB001EN in case you need to remove and re-mount the syringes of the titration unit.

6.9 Venting the Gas Supply Line (not applicable to SurPASS 3 Eco)

Proceed as follows to disconnect the SurPASS 3 instrument from your in-house gas supply:

1. Make sure the external gas supply is activated in the software.
2. Switch off the nitrogen supply in your laboratory.
3. Watch the display on the manometer of your in-house gas supply and wait for the pressure to drop. This might take 1-2 minutes.
4. Switch off the instrument by using the power switch on the rear side of the SurPASS 3 instrument.
5. Detach the pressure hose (hose (5) in Fig. 18) by pressing down and holding its releasing ring. Remove the hose.

6.10 Storing and Transporting the Instrument

- Store the instrument under clean conditions that fulfill the ambient conditions provided in appendix A Technical Data.

- Make sure that it was rinsed with ultra-pure water and that it is free of salt and aggressive chemical liquids.
- All liquids and samples residues must be removed from the measuring cells if not in use.
- Empty and rinse the acid and base titration lines to keep up the titration syringes if the instrument is not in use.
- Turn off the in-house nitrogen supply and vent the gas supply line. Refer to section 6.9 Venting the Gas Supply Line (not applicable to SurPASS 3 Eco) for details.
- Empty the measuring cell before you move or lift the instrument.
- Store the instrument in a dry and empty condition.
- Store the pH electrode in the appropriate storage beaker filled with 3 mol/L KCl.
- Store the conductivity electrode in the storage beaker of the instrument in dry condition.
- The SurPASS 3 must be transported in the original package or in the corresponding transport box.
- During transport, the instrument must not be exposed to intense shocks or high forces.

6.11 Packing the Instrument for Returns

This should be performed by a qualified service technician. Contact your local Anton Paar representative for details.

TIP: Find the contact data of your local Anton Paar representative on the Anton Paar website (<http://www.anton-paar.com>) under "Contact".

7 Maintenance and Repair

7.1 Maintenance Performed by an Authorized Anton Paar Service Engineer

The SurPASS 3 requires a periodical maintenance which shall be performed by an authorized Anton Paar Service Engineer¹.

A missing maintenance may mean that under certain conditions your warranty is no longer valid².

Maintenance Interval:

- Once a year,
- Or if the standard maintenance does not lead to a successful verification,
- Whichever is sooner.

Parts to be exchanged at every service interval (wear and tear parts):

- Set of O-rings for the plug-in module (including hydraulic valve block, pressure vessel, front panel and hydraulic valve motor)
- O-rings of the electrodes
- Titration hoses and titration syringes if corroded
- Electrolyte hoses
- Nitrogen purge tube and throttle
- Gaskets and O-rings of the measuring cells

Parts to be controlled at every service interval and to be exchanged if required (wear and tear parts):

- Ceramic valve disks
- Membrane pump
- Digital pressure regulator
- Conductivity electrode
- pH electrode

7.2 Repair Performed by an Authorized Anton Paar Representative

In case your instrument needs repair, contact your local Anton Paar representative, who will take care of the necessary steps.

If your instrument needs to be returned, request an RMA (Return Material Authorization Number). It must not be sent without the RMA and the filled "Safety Declaration for Instrument Repairs". Please make sure it is cleaned before return.

TIP: Find the contact data of your local Anton Paar representative on the Anton Paar website (<http://www.anton-paar.com>) under "Contact".

¹ Please contact your Anton Paar representative to get an offer.

² For detailed information please see the general terms of delivery (GTD) on the Anton Paar website (<http://www.anton-paar.com>).

Appendix A: Technical Data

Measuring range

Streaming potential	$\pm 2000 \text{ mV} \pm (0.2 \% + 4 \text{ } \mu\text{V})$
Streaming current	$\pm 2 \text{ mA} \pm (0.2 \% + 1 \text{ pA})$
Cell resistance	$5 \text{ } \Omega \dots 20 \text{ M}\Omega \pm (2 \% + 0.5 \text{ } \Omega)$
Differential pressure	SurPASS 3: 3500 mbar $\pm (0.2 \% + 0.5 \text{ mbar})$ external pressure supply (compressed N ₂) required for differential pressure > 1500 mbar. SurPASS 3 Eco: 1200 mbar $\pm (0.2 \% + 0.5 \text{ mbar})$
Conductivity	$0.1 \dots 2 \times 10^5 \text{ mS/m}$
pH measurement	pH 1 ... 14 ± 0.05
Temperature	$5^\circ \dots 40^\circ \text{ C} \pm 1.5^\circ \text{ C}$

Electrolyte solution (reasonable measuring range for Ag/AgCl electrodes)

pH value	pH 2 ... 12
Conductivity	0.1 ... 1000 mS/m
Temperature	20 ... 40 °C
Water	ASTM I grade

Electrolyte solution (reasonable measuring range for gold electrodes)

pH value	pH 2 ... 14
Conductivity	of 0.1 ... 150 mS/m for fiber, powder and conductive samples
Conductivity	of 0.1 ... 700 mS/m for planar, inert and not conductive samples
Temperature	20 ... 40 °C
Water	ASTM I grade

Mains supply

Power consumption	200 VA
Fuses	T4A H 250V

Dimensions

Footprint	$612 \times 536 \text{ mm (D} \times \text{W)}$
-----------	---

Appendix A: Technical Data

Clearance	360 mm
Weight	SurPASS 3: 26 kg SurPASS 3 Eco: 24 kg

Environmental conditions

Ambient temperature	20 ... 40 °C
Humidity	10 ... 80 % r.h. not condensing
Overvoltage category	II according to EN 61010-1
Pollution degree	2 according to EN 61010-1
Noise emission	30 dB

Sample size requirement

Clamping Cell	min. 35 mm × 15 mm or min. 17 mm in diameter, max. thickness 40 mm
Cylindrical Cell	particle size > 25 µm
Adjustable Gap Cell	20 mm × 10 mm, max. thickness 2 mm
Adjustable Gap Cell for Disks	14 mm or 15 mm in diameter, max. thickness 2 mm
Measuring Cell for Ceramic Membrane 25 mm	Diameter: 25.2 ± 0.3 mm; Length: 100 mm ± 4 mm Diameter: 10 ± 0.1 mm; Length: 96 mm ± 4 mm
Measuring Cell for Ceramic Membrane 30 mm	Diameter: 30.8 ± 0.3 mm; Length: 100 mm ± 4 mm Diameter: 13.1 ± 0.1 mm; Length: 96 mm ± 4 mm

Computer specification

Operating system	Microsoft Windows® 7 32 bit and 64 bit Service Pack 1 or later (Windows 7 Starter Edition is not supported) Microsoft Windows® 8 32 bit and 64 bit Microsoft Windows® Server 2008 / 2008 R2 32 bit and 64 bit Service Pack 2 or higher for Windows Server 2008 Microsoft Windows® 10 32 bit and 64 bit
CPU	Intel® Core™ i3
RAM	4 GB (depends on operating system)
Hard disk	20 GB free disk space
Video	Screen resolution at least 1024 x 768 at True Color (32 bit)
Communication port	1 free USB port

The SurPASS 3 is an instrument for indoor use at an altitude of up to 2500 m.

Appendix B: Declaration of Conformity

EC Declaration of Conformity



The manufacturer Anton Paar GmbH, Anton-Paar-Str. 20, A-8054 Graz, Austria – Europe, hereby declares that the machinery described below:

Description: **SURPASS 3 SURFACE POTENTIAL ANALYZER FOR SOLID SAMPLES**
 Model: **SURPASS 3**
 Material number: **20880, 45030**
 Serial number:

complies with all the relevant provisions of the Machinery Directive (2006/42/EC, O.J. L 157/24 of 9.6.2006) and the regulations transposing it into national law

complies with all the relevant provisions of the Electromagnetic Compatibility Directive (2014/30/EU, O.J. L 96/79 of 29.3.2014)

complies with all the relevant provisions of the RoHS Directive (2011/65/EU, O.J. L 17/488 of 1.7.2011)

complies with all the safety objectives of the Low Voltage Directive (2014/35/EU, O.J. L 96/357 of 29.3.2014)

complies with the provisions of the following harmonized standards:

- EN ISO 12100:2010 Safety of machinery - General principles for design - Risk assessment and risk reduction
 - EN 61326-1:2013 Electrical equipment for measurement, control and laboratory use - EMC requirements - Part 1: General requirements
- complies with the provisions of the following technical standards:
- EN 61010-1:2010 +A1:2019 Safety requirements for electrical equipment for measurement, control and laboratory use – Part 1: General requirements

This declaration relates exclusively to the machinery in the state in which it was placed on the market, and excludes components which are added and/or operations carried out subsequently by the final user.

The manufacturer complies the technical file according to 2006/42/EC Annex II

Place and date of issue: Graz, 2022-08-01


 DI Alfred Freiberger
 Executive Director
 Business Unit Characterization


 DI Dr. Andreas Kautsch
 Head of Particle Analysis & Surface Charge
 Business Unit Characterization

UK Declaration of Conformity



The Manufacturer Anton Paar GmbH, Anton-Paar-Str. 20, A-8054 Graz, Austria – Europe hereby declares that the product listed below

Description: **SURPASS 3 SURFACE POTENTIAL ANALYZER FOR SOLID SAMPLES**
 Model: **SURPASS 3**
 Material number: **20880, 45030**
 Serial number:

is in conformity with all the relevant UK legislation

Supply of Machinery (Safety) Regulations 2008, 2008 No. 1597

Electromagnetic Compatibility Regulations 2016, 2016 No. 1091

Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic Equipment Regulations 2012, 2012 No. 3032

is in conformity with all the UK safety objectives of the

Electrical Equipment (Safety) Regulations 2016, 2016 No. 1101

complies with the designated standards:

- EN ISO 12100:2010
- EN 61326-1:2013
- EN 61010-1:2010 +A1:2019 +A1:2019/AC:2019

The declaration of conformity is issued under the sole responsibility of the manufacturer.

Importer: Anton Paar Ltd, Unit F, The Courtyard, Hatfield Rd, St Albans AL4 0LA, United Kingdom.

Place and date of issue: Graz, 2022-08-01


 DI Alfred Freiberger
 Executive Director
 Business Unit Characterization


 DI Dr. Andreas Kautsch
 Head of Particle Analysis & Surface Charge
 Business Unit Characterization

Appendix C: Requirements for Nitrogen Supply



WARNING

Risk of injury

As there is a risk of suffocation when using nitrogen for measurement, ensure adequate ventilation of the room.

External nitrogen supply is required for high pressure application (> 1500 mbar), adsorption studies or purging the electrolyte solution to prevent carbon dioxide dissolution.

The following requirements need to be fulfilled:

- Use nitrogen 5.0 for any of these purposes.
- If nitrogen is supplied directly from a gas container, a double-stage gas valve is required to reduce the primary pressure.
- If nitrogen is supplied from an in-house gas supply, where the primary pressure is already reduced, a single-stage gas valve is required.
- The operating pressure, i.e. the pressure applied on the position "IN" of the pressure control unit, shall be set to 6 bar to ensure correct functioning of the pressure control unit.
- The maximum operating pressure must not exceed 10 bar.
- The outlet port of the external gas supply should accommodate a pressure hose with 6 mm (1/4 inch).

NOTICE

The maximum pressure applicable to the instrument is 3500 mbar (high-pressure application).



Fig. C-1: Air Liquide pressure-reducing valve BS 20-1-2 with a resolution of 0.05 bar for use with a nitrogen supply with 20 bar primary pressure.

NOTICE

If using an inert gas instead of Nitrogen e.g. argon, make sure that the quality of the gas is 5.0.

The gas consumption of SurPASS 3 during operation is as follows:

- 0.2 l/min for purging the electrolyte solution
- 0.3 l/min for rinse and measurement cycles.